

Quick-Start Guide MIKRO 2.0 | 2.0R


This Quick-Start Guide only serves as a reference for the initial set up of the machine covering the most important functions and operating steps. It does not replace the full operating instructions. Please read the manual, including all safety instructions before use.

Initial Setup


1. Switch on the device.




2. Remove the transport protection from the rotor.

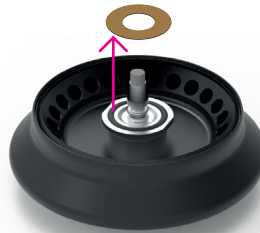


3. Place the rotor lid vertically onto the rotor and push down firmly. The removal aid and unlocking pin must be flush.

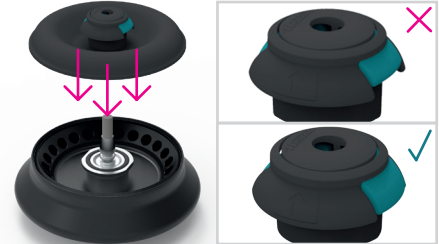

4. To open the lid of the centrifuge press

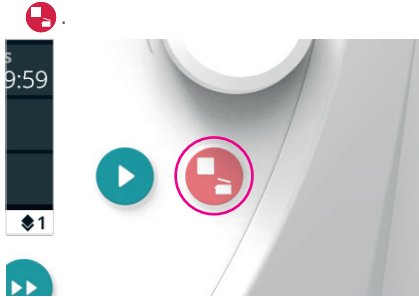


5. Install the rotor vertically.

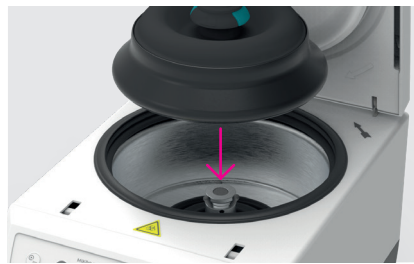

6. To close the lid, press the front edge of the lid down slightly.



7. Functions of the TURN/PUSH button. Select and confirm.

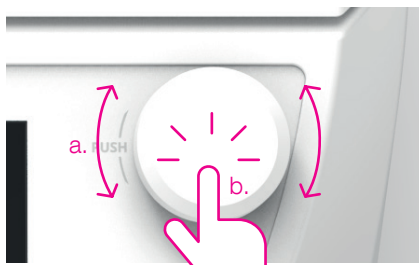


7. a. Turn = adjust parameters.




7. b. Press = confirm parameter value and/or navigate to the next parameter.



8. a. To open the sub menu press . b. To select P5 to P8, hold and select your program button.

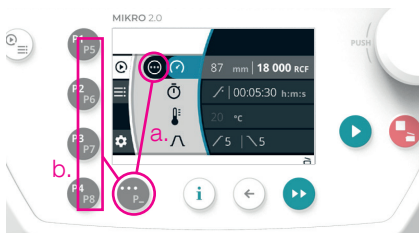

9. To switch between the Quick-Adjust and Program windows press . To close notifications and go back, press .

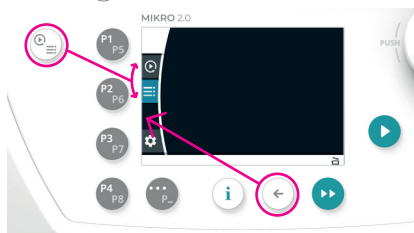


10. Press to start centrifugation.




Remove Rotor


1. To stop, press . For quick stop press the button twice. To open the lid press again.

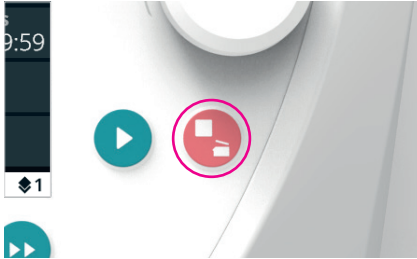

2. a. Press the release pin down with your thumb. b. reach under the knob with your index and middle finger. Pull the rotor out.

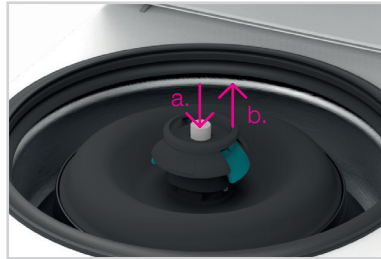

3. Press the lid release together with your index finger and thumb and lift the lid.

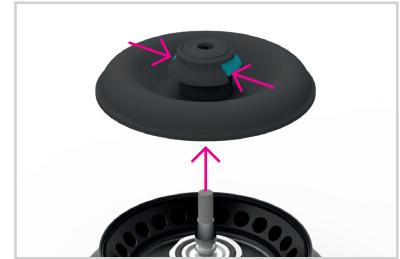


Create and start an individual program (authorization level 3)


1. Press to switch to Program window.




2. Navigate with the TURN/PUSH-button to .

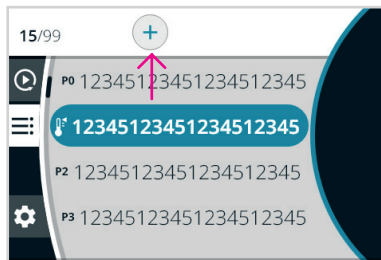


3. Press the TURN/PUSH-button to select „add new program“.

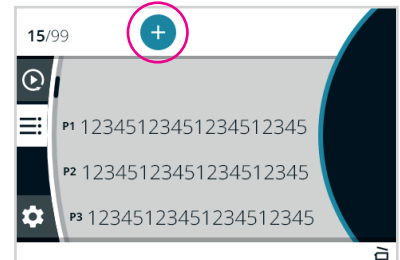


4. Create a program name using the TURN/PUSH button to navigate the keypad and select letters.

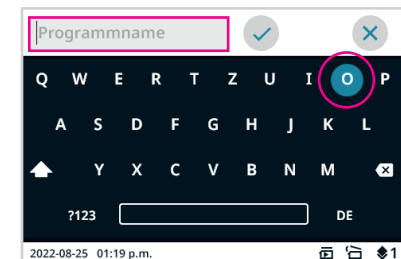


5. Press after entering the program name to leave the keypad.

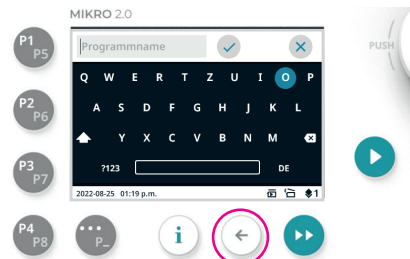


6. Save the program name by pressing the TURN/PUSH button. To discard, turn and then press the button.

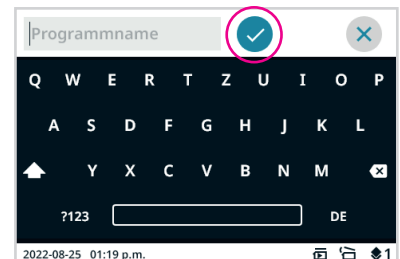

7. Select the new program by press the TURN/PUSH button. Press the TURN/PUSH again to adjust the parameters.

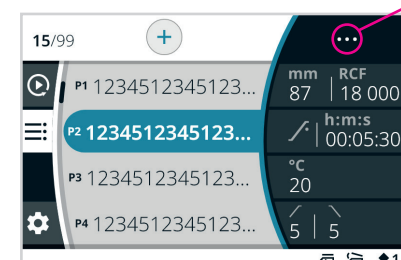

8. Press to open the program sub menu. (a.) Set as a favorite (b.) Copy (c.) Change name (d.) Delete

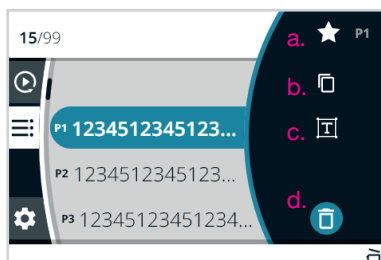

9. Press to return to the program window. To start the program press .

