EBA 270

Inhalt des Dokuments / content of the document

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Operating instructions (EN)

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Operating instructions

EBA 270

Translation of the original operating instructions

©2023 – All rights reserved

Andreas Hettich GmbH & Co. KG

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D-78532 Tuttlingen, Germany

Telephone: +49 (0)7461 705-0

Fax: +49 (0)7461 705-1125

Email: info@hettichlab.com, service@hettichlab.com

Internet: www.hettichlab.com

Table of contents

| | | |
|----------|---|-----------|
| 1 | About this document. | 5 |
| 1.1 | Use of this document. | 5 |
| 1.2 | Gender reference. | 5 |
| 1.3 | Symbols and labels in this document. | 5 |
| 2 | Safety. | 5 |
| 2.1 | Intended use. | 5 |
| 2.2 | Personnel requirements. | 6 |
| 2.3 | Operator's responsibility. | 7 |
| 2.4 | Safety instructions. | 7 |
| 3 | Device overview. | 9 |
| 3.1 | Technical data. | 9 |
| 3.2 | European registration. | 11 |
| 3.3 | Important labels on the packaging. | 11 |
| 3.4 | Important labels on the device. | 12 |
| 3.5 | Operating and indicator elements. | 13 |
| 3.5.1 | Control. | 13 |
| 3.5.2 | Indicator elements. | 13 |
| 3.5.3 | Controls. | 13 |
| 3.6 | Original spare parts. | 14 |
| 3.7 | Scope of delivery. | 14 |
| 3.8 | Returns. | 14 |
| 4 | Transport and storage. | 15 |
| 4.1 | Transport and storage conditions. | 15 |
| 4.2 | Fastening the transport lock. | 15 |
| 5 | Commissioning. | 16 |
| 5.1 | Unpacking the centrifuge. | 16 |
| 5.2 | Removing the transport lock. | 17 |
| 5.3 | Setting up and connecting the centrifuge. | 17 |
| 5.4 | Switching the centrifuge on and off. | 19 |
| 6 | Operation | 19 |
| 6.1 | Opening and closing the lid. | 19 |
| 6.2 | Removing and installing the rotor. | 19 |
| 6.3 | Loading. | 20 |
| 6.4 | Centrifugation. | 21 |
| 6.4.1 | Centrifugation in continuous operation. | 21 |
| 6.4.2 | Centrifugation with time preselection. | 22 |
| 6.4.3 | Short-time centrifugation. | 22 |
| 7 | Software operation. | 23 |
| 7.1 | Centrifugation parameters. | 23 |
| 7.1.1 | Set the brake level. | 23 |
| 7.1.2 | Relative centrifugal force, RCF. | 23 |

| | | |
|-----------|--|-----------|
| 7.1.3 | Centrifugation of substances or mixtures of substances with a density higher than 1.2 kg/dm ³ . | 23 |
| 7.1.4 | Centrifuging radius. | 24 |
| 7.2 | Machine Menu. | 24 |
| 7.2.1 | Audible signal. | 24 |
| 7.2.1.1 | General. | 24 |
| 7.2.1.2 | Setting an audible signal. | 24 |
| 8 | Cleaning and care. | 25 |
| 8.1 | Overview table. | 25 |
| 8.2 | Instructions for cleaning and disinfection. | 25 |
| 8.3 | Cleaning. | 26 |
| 8.4 | Disinfection. | 26 |
| 8.5 | Maintenance. | 27 |
| 9 | Troubleshooting. | 28 |
| 9.1 | Fault description. | 28 |
| 9.2 | Performing a MAINS RESET. | 29 |
| 9.3 | Emergency release. | 29 |
| 9.4 | Replacing the mains input fuse. | 30 |
| 10 | Disposal. | 31 |
| 10.1 | General instructions. | 31 |
| 11 | Index. | 32 |

1 About this document

1.1 Use of this document

- Read this document carefully and in full before commencing initial operation of the device.
Observe other enclosed instruction sheets where necessary.
- This document constitutes part of the device and must be kept within easy reach.
- This document must be included if the device is passed on to a third party.
- The most recently updated version of this document in the available languages can be found on the manufacturer's website at: ➔ <https://www.hettichlab.com/de/download-center/>

1.2 Gender reference

The employed masculine or feminine language form is to facilitate reading. In the spirit of equal treatment, corresponding terms apply in principle to all genders and do not imply any valuation.

1.3 Symbols and labels in this document

General symbols

The following markers are used in this document to highlight instructions, results, listings, references and other elements:

| Marker | Explanation |
|---|---|
| 1.  2.  3.  ...  | Step-by-step instructions |
|  | Results of action steps |
|  | References to sections of the document and other applicable documents |
| ■ ... ■ ... | Listings without a fixed order |
| <i>[Buttons]</i> | Controls (for example: buttons, switches) |
| <i>'Indicator'</i> | Indicator elements (for example: signal lights, screen elements) |

2 Safety

2.1 Intended use

Intended use

The centrifuge **EBA 270** is an in vitro diagnostic medical device in accordance with the In Vitro Diagnostic Medical Devices Regulation (EU) 2017/746. The device is used for centrifugation as well as enrichment of sample material of human origin for subsequent further processing for diagnostic purposes. The user can set each of the variable physical parameters within the limits set by the device.

The centrifuge may only be used by qualified personnel in closed laboratories. The centrifuge is only intended for the use referred to above. Intended use also includes observing all instructions in the user manual and compliance with inspection and maintenance. Any other use or use beyond this is considered improper. Andreas Hettich GmbH & Co. KG shall not be liable for any damage arising from this.

Non-intended use

- The centrifuge is not suitable for use in explosive or radioactive, or biologically or chemically-contaminated atmospheres.
- The user must take appropriate actions when centrifuging hazardous substances or mixtures of substances that are toxic, radioactive or contaminated with pathogenic microorganisms.
The manufacturer generally recommends using only centrifuge tubes with special screw caps designed for use with hazardous substances. Use sealable centrifuge tubes with a biosafety system for materials of risk groups 3 and 4.
- The manufacturer does not recommend centrifugation of flammable or explosive materials.
- The manufacturer does not recommend centrifugation of materials that react chemically with one another with high activation energy.

Foreseeable misuse

The manufacturer recommends using only accessories approved that it has approved for the intended purpose.

Only operate the centrifuge under supervision.

2.2 Personnel requirements

Required qualifications

The user has read the Operating Manual in full and familiarised themselves with the device.

NOTICE

Damage to the device by unauthorised personnel

- Tampering with and modifications to devices by unauthorised persons are at the operating organisation's own risk and will result in the loss of all warranty and liability claims.

Trained user

The user has been educated and trained in laboratory work and is able to carry out the work assigned to them, and to recognise and prevent potential hazards independently.

Personal protective equipment

Lack of personal protective equipment or unsuitable personal protective equipment increases the risk of impaired health and injury.

- Only use personal protective equipment that is in proper condition.
- Only use personal protective equipment that is adapted to the person (correct size, for example).
- Observe instructions on other protective equipment for specific activities.

2.3 Operator's responsibility

Follow the instructions in this document for proper and safe use of the device.

Keep the user manual for future reference.

Provide information

- Following the instructions in this document will help:
 - To avoid dangerous situations.
 - To minimise repair costs and downtime.
 - To increase the reliability and service life of the device.
- The operator is responsible for compliance with company regulations, standards and national laws.
- Note and keep the revision of the document separate from the document. If lost, the document can be replaced in the correct revision.
- Keep the user manual available at the place where the device is used.
- Pass the user manual on to the buyer when the device is sold.

Personnel training

Lack of knowledge when working with the device may result in serious injury or death.

- Instruct personnel on their tasks and the associated risks in accordance with the instruction.

2.4 Safety instructions

Reporting serious incidents and notifiable incidents

In the event of serious incidents or notifiable incidents involving the device or its accessories, these must be reported to the manufacturer and, where applicable, to the competent authority where the user and/or the patient is registered.

DANGER

Risk of contamination for the user due to inadequate cleaning or failure to observe the cleaning instructions.

- Observe cleaning instructions.
- Wear personal protective equipment when cleaning the device.
- Observe laboratory regulations (e.g. TRBAs, the German Protection against Infection Act, hygiene plan) for handling biological agents.

DANGER

Fire and explosion hazard due to hazardous substances in samples.

- Observe relevant regulations and directives for handling chemicals and hazardous substances.
- Do not use aggressive chemicals (for example: dangerous, corrosive extraction agents such as chloroform, strong acids).

**WARNING**

Dangers due to insufficient maintenance or maintenance not carried out on time.

- Follow maintenance intervals.
- Check the device for visible damage or defects.
If any visible damage or defects are present, take the device out of service and inform a service technician.

 **WARNING**

Risk of electric shock due to ingress of water or other liquids.

- Protect the device against external liquids.
- Do not pour any liquids into the interior of the device.
- Transport using original transport packaging.

 **WARNING**

Contamination with hazardous substances and substance mixtures!

Observe the following actions for substances and substance mixtures that are toxic, radioactive and/or contaminated with pathogenic microorganisms:

- As a rule, use only centrifuge tubes with special screw caps for hazardous substances.
- Use sealable centrifuge tubes with a biosafety system for materials of risk groups 3 and 4.
- If no biosafety system is used, the device is not micro-biologically tight in the sense of standard EN / IEC 61010-2-020.
- Contact the manufacturer if necessary.

**WARNING**

Risk of injury and damage to the device due to a loose rotor.

- The driver of the rotor shaft must be correctly seated in the groove of the rotor when mounting the rotor.
- Hand-tighten the nut securing the rotor.
- Check that the rotor is firmly seated.
- Follow maintenance intervals.

**CAUTION**

Risk of injury due to rotating rotor

Long hair and items of clothing can get caught on the rotor if the rotor is moved manually.

- Tie long hair back.
- Do not allow garments to hang in the centrifuging chamber.

NOTICE

Damage to the device electronics due to incorrect voltage or frequency at the device circuit breaker.

- Operate the device with the correct mains voltage and mains frequency.
The value can be found in the technical data and on the rating plate.

NOTICE

Damage to the device and samples due to premature program termination.

Premature program termination is caused by power failure, switching off during the program or pulling out the mains plug.

- Do not switch off the device while the program is running.
- Do not trigger the emergency release on the device while the program is running.
- Do not pull out the mains plug while the program is running.

3 Device overview

3.1 Technical data

| | | |
|---|--|--------------|
| Manufacturer | Andreas Hettich GmbH & Co. KG, D-78532 Tuttlingen | |
| Model | EBA 270 | |
| Type | 2300 | 2300-01 |
| Mains voltage ($\pm 10\%$) | 200-240 V 1~ | 100-127 V 1~ |
| Mains frequency | 50-60 Hz | 50-60 Hz |
| power consumption | 130 VA | 125 VA |
| Power consumption | 0.7 A | 1.25 A |
| max. capacity | 16 x 15 ml | |
| max. permissible density | 1.2 kg/dm ³ | |
| max. speed (RPM) | 4000 | |
| max. acceleration (RCF) | 2254 | |
| max. kinetic energy | 250 Nm | |
| Obligation to perform checks (DGUV Rules 100-500) (valid only in Germany) | No | |

| | | |
|--|---|-------------|
| Ambient conditions (EN / IEC 61010-1): | | |
| Installation site | indoors only | |
| Altitude | up to 2000 m above sea level | |
| Ambient temperature | 2 °C to 40 °C | |
| Humidity | maximum relative humidity 80% for temperatures up to 31 °C, decreasing linearly to 50% relative humidity at 40 °C. | |
| Overvoltage category (IEC 60364-4-443) | II | |
| Pollution level | 2 | |
| Device protection class | I not suitable for use in potentially explosive atmospheres. | |
| EMC: | | |
| Emitted EM interference, EM interference immunity | EN / IEC 61326-1 Class B | FCC Class B |
| Noise level (rotor-dependent) | ≤51 dB(A) | |
| Dimensions: | | |
| Width | 326 mm | |
| Depth | 389 mm | |
| Altitude | 239 mm | |
| Weight | approx. 13.5 kg | |

Rating plate

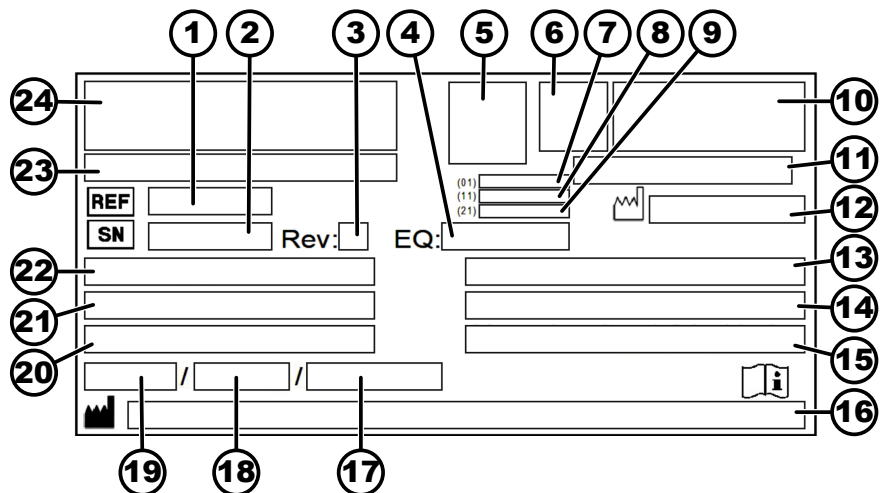
Fig. 1: Rating plate

- 1 Item number
- 2 Serial number
- 3 Revision
- 4 Equipment number

- 5 Data matrix code
- 6 any labelling indicating whether medical device or in vitro diagnostic medical device
- 7 Global Trade Item Number (GTIN)
- 8 Date of manufacture
- 9 Serial number
- 10 any EAC mark, CE mark
- 11 Country of manufacture
- 12 Date of manufacture
- 13 Mains frequency
- 14 Maximum kinetic energy
- 15 Maximum permissible density
- 16 Manufacturer's address
- 17 any Coolant circuit pressure
- 18 any Coolant capacity
- 19 any Coolant type
- 20 Revs per minute
- 21 Performance values
- 22 Mains voltage
- 23 any Device designation
- 24 Manufacturer's logo

3.2 European registration

Device conformity

Device conformity according to EU directives.

Single Registration Number

SRN: DE-MF-000010680

Basic-UDI-DI

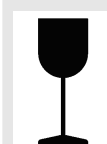
| Basic-UDI-DI | Device assignment |
|------------------|--|
| 040506740100079W | EBA 270 (in vitro diagnostic medical device) |

3.3 Important labels on the packaging

TOP

This is the correct upright position of the shipping container for transport and/or storage.

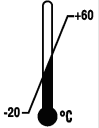
FRAGILE GOODS

The contents of the shipping container are fragile, so it must be handled with care.

PROTECT FROM MOISTURE
The shipping container must not be exposed to rain and kept in a dry environment.

TEMPERATURE LIMITATION
The shipping container must be stored, transported and handled within the indicated temperature range (-20 °C to +60 °C).

HUMIDITY LIMITATION
The shipping container must be stored, transported and handled within the indicated humidity range (10% to 80%).

STACK LIMITATION BASED ON QUANTITY
Maximum number of identical packages that may be stacked on the lowest package, "n" standing for the number of packages allowed. The lowest package is not included in "n".

3.4 Important labels on the device

The signs on the device must not be removed or covered, or have anything pasted over them.

Attention, general danger area.
Ensure you read the instructions for commissioning and operation and observe the safety instructions before using the device.

Biohazard warning.

Direction of rotation of the rotor.
The orientation of the arrow indicates the rotor's direction of rotation.

Direction of rotation of the emergency release.

Symbol for the separate collection of electrical and electronic equipment, in accordance with Directive 2012/19/EU (WEEE).

Use in European Union countries, Norway and Switzerland.

3.5 Operating and indicator elements

3.5.1 Control

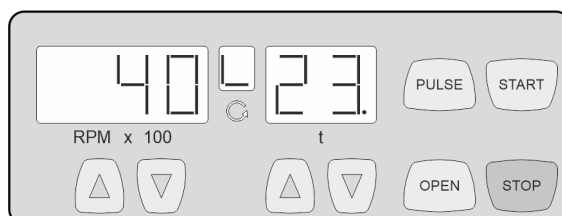
Fig. 2: Control

3.5.2 Indicator elements

Fig. 3: 'Lid closed' indicator

Fig. 4: 'Lid open' indicator

Fig. 5: 'Rotation' indicator

3.5.3 Controls

Fig. 6: [Mains switch]

Fig. 7: [OPEN] button

Fig. 8: [PULSE] button

Fig. 9: [START] button

- The indicator appears when the lid is closed.
- If the 'Lid closed' and 'Lid open' indicators flash alternately, further operation of the centrifuge is only possible after opening the lid once.
- The indicator appears when the lid is open.
- The indicator light rotates when the rotor is turning.
- Switch the device on and off.
- Unlock the lid.
- Short-time centrifugation.
- The centrifugation run takes place as long as the button is being pressed.
- Indicator brake level.
- Start centrifugation run.

Fig. 10: [STOP] button

RPM x 100

Fig. 11: [RPMx100] buttons

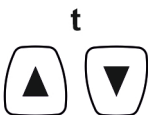
Fig. 12: [t] buttons

- End the centrifugation run.
The rotor ramps down to a stop at the preselected brake level.
- Save brake level.
- A numerical value from 500 RPM to the maximum rotor speed can be set.
Adjustable in increments of 100 (RPM = displayed value x 100).
- Indicator brake level.
- Enter runtime.
- Brake level 0 or 1.

3.6 Original spare parts

Use only original spare parts from the manufacturer and approved accessories.

3.7 Scope of delivery

The following accessories are supplied with the centrifuge:

- 2 Fuse link
- 1 Release pin
- 1 single-end open-ended spanner (10 mm AF)
- 1 right-angled hex key (5 mm)
- 1 hex key (SW5 x 100)
- 1 power cable
- 1 Rotor
- 6 sleeves, 10 ml
- 6 sleeves, 15 ml
- 1 user manual
- 1 instruction sheet, transport lock

Rotors and the corresponding accessories are supplied depending on the order.

3.8 Returns

An original Return Material Authorisation (RMA) form from the manufacturer must always be requested for a return. Secure and reliable acceptance and booking in of the goods with the manufacturer is not possible without an original RMA form from the manufacturer. The Return Material Authorisation (RMA) form contains a Declaration of No Objection (UBE), which must be completed in full and enclosed with the return.

If the device and/or accessories are returned to the manufacturer, the complete return shipment must be cleaned and decontaminated by the sender. If returns are not cleaned and/or decontaminated or are insufficiently cleaned and/or decontaminated, this will be performed by the manufacturer and charged to the sender.

The original transport locks must be attached for return shipment, see [↗ Chapter 4 'Transport and storage' on page 15](#). The device must be shipped in its original packaging.

4 Transport and storage

4.1 Transport and storage conditions

Transport conditions

NOTICE

Damage to the device due to failure to use the transport locks.

- Secure the transport locks before transporting the device.

NOTICE

Damage to the device due to condensation.

There is a risk of condensation forming on electrical components when component surfaces are cold and the surrounding air is warmer. The condensation that forms may cause a short circuit and/or destroy electronics.

- Warm the device up for at least 3 hours in a warm room before connecting it to the mains.
or
- Warm up for 30 minutes in a cold room.

- Before transporting, fasten the transport lock and disconnect the device from the mains socket.
- The transport temperature must be between -20 °C and +60 °C.
- Humidity must not be condensing. Humidity must be between 10% and 80%.
- Be aware of the weight of the device.
- When transporting using a transport aid (e.g., a pallet truck), the transport aid must be able to carry at least 1.6 times the transport weight of the device.
- Secure the device to prevent it tipping over and falling down during transport.
- Never transport the device sideways or upside down.

Storage conditions

- The device must be stored in the original packaging.
- Only store the device in dry rooms.
- The storage temperature must be between -20 °C and +60 °C.
- Humidity must not be condensing. Humidity must be between 10% and 80%.

4.2 Fastening the transport lock

Personnel:

- Trained user

The lid is closed.
 The mains cable is disconnected from the device.

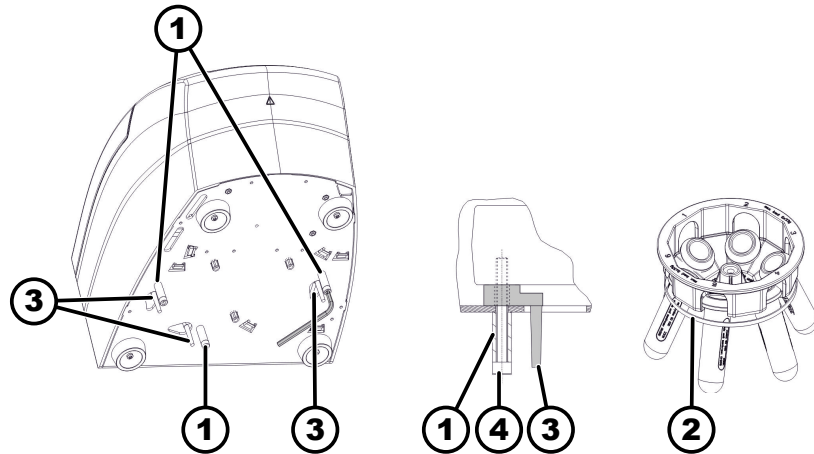
Fig. 13: Transport lock

- 1 Spacer sleeves
- 2 Rubber band
- 3 Transport lock
- 4 Screws

1. ➤ Tilt the device on the back of the device.
2. ➤ Insert 3 transport locks (3).
3. ➤ Screw in 3 screws (4) with spacer sleeves (1).
4. ➤ Secure the rotor bucket using a rubber band (2).

5 Commissioning

5.1 Unpacking the centrifuge

CAUTION

Danger of crushing due to parts falling out of the transport packaging.

- Keep the device balanced during the unpacking process.
- Only open the packaging at the points provided for this purpose.

CAUTION

Risk of injury from lifting heavy loads.

- Provide an adequate number of helpers.
- Note the weight. See ➔ Chapter 3.1 'Technical data' on page 9.

NOTICE

Damage to the device due to improper lifting.

- Do not lift the centrifuge by the control panel or the control panel holder.

Personnel:

- Trained user

1. ➤ Open the box at the top.
2. ➤ Remove the padding.
3. ➤ Remove the device and accessories by lifting them up out of the box.
4. ➤ Place the device on a stable and level surface.

5.2 Removing the transport lock

Personnel:

- Trained user

The lid is closed.

The mains cable is disconnected from the device.

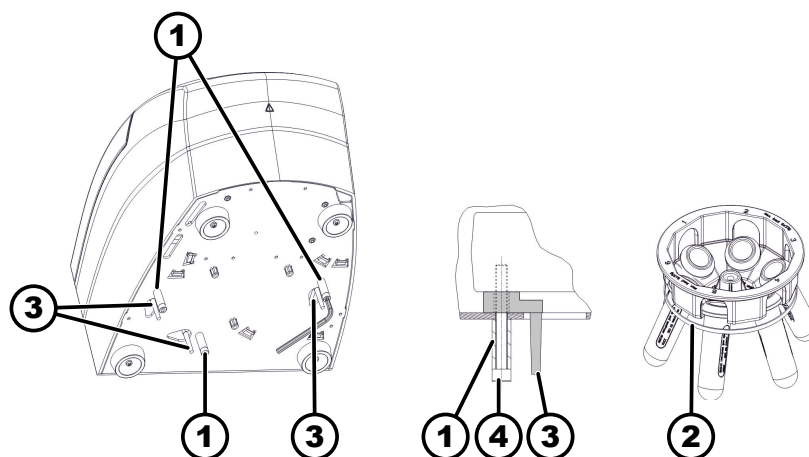
Fig. 14: Transport lock

- 1 Spacer sleeves
- 2 Rubber band
- 3 Transport lock
- 4 Screws

1. ➤ Tilt the device on the back of the device.
2. ➤ Unscrew 3 screws (4) with spacer sleeves (1).
3. ➤ Remove 3 transport locks (3).
4. ➤ Keep the screws, spacer sleeves and transport lock in a safe place.
5. ➤ Remove the rubber band (2) from the rotor.

5.3 Setting up and connecting the centrifuge

Setting up the centrifuge

WARNING

Risk of injury due to failing to maintain a sufficient distance to the centrifuge.

- As per EN / IEC 61010-2-020, no persons, hazardous materials or objects may be present within a **safety zone of 300 mm** around the centrifuge during a centrifugation run.
- A distance of **300 mm** from the ventilation slots and ventilation openings of the centrifuge must be maintained.

**CAUTION**

Risk of crushing and damage to the device due to it falling down because of vibration-induced position alterations.

- Place the device on a stable and level surface.
- Select the installation surface dependent on the weight of the device.

**NOTICE**

Damage to the samples and the device if the ambient temperature exceeds or falls below the respective maximum/minimum permissible ambient temperature.

- Comply with the maximum and minimum permissible ambient temperatures for installation of the device.
- Do not place the device next to a heat source.
- Do not expose the device to direct sunlight.
- Do not expose the device to frost.

Personnel:

- Trained user

1. Place the device on a stable and level surface.
2. Maintain a distance of 300 mm around the device.
3. Comply with the ambient conditions in the technical data (→ *Chapter 3.1 'Technical data' on page 9*).

Connecting the centrifuge**NOTICE**

Damage to the device by unauthorised personnel

- Tampering with and modifications to devices by unauthorised persons are at the operating organisation's own risk and will result in the loss of all warranty and liability claims.

**NOTICE**

Damage to the device due to condensation.

There is a risk of condensation forming on electrical components when component surfaces are cold and the surrounding air is warmer. The condensation that forms may cause a short circuit and/or destroy electronics.

- Warm the device up for at least 3 hours in a warm room before connecting it to the mains.
or
- Warm up for 30 minutes in a cold room.

Personnel:

- Trained user

1. A type B residual current circuit breaker must be used if the device is additionally protected with a residual current circuit breaker in the building installation.

When using a different type, the residual current circuit breaker may either not switch off the unit if there is a fault on the unit, or it may switch off the unit even though there is no fault on the unit.

2.  Check whether the mains voltage matches the specification on the rating plate.
3.  Connect the device to a standard mains socket using the mains cable.

5.4 Switching the centrifuge on and off.

Switching the centrifuge on

Personnel:

- Trained user
-  Set the mains switch to *///*.
- The buttons flash, depending on the centrifuge type.
The last centrifugation data used is displayed.

Switching off the centrifuge

The rotor is stationary.

-  Set the mains switch to */0/*.

6 Operation

6.1 Opening and closing the lid

Opening the lid

Personnel:

- Trained user
- The centrifuge is switched on.
The rotor is stationary.
1.  Press the *[OPEN]* button.
 - The 'Lid open' indicator appears.
 2.  Open the lid.

Closing the lid

NOTICE

Damage to the device caused by the lid slamming.

- Close the lid slowly.
- Do not slam the lid.

Personnel:

- Trained user
-  Close the lid and press the front edge of the lid down gently.
 - The 'Lid closed' indicator appears.

6.2 Removing and installing the rotor

Removing the rotor

Personnel:

- Trained user

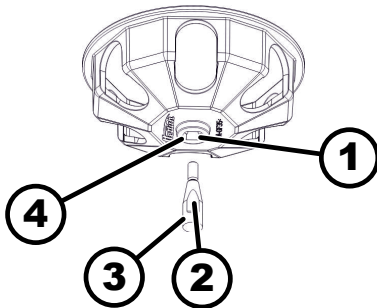
Fig. 15: Rotor installation and removal

- 1 Groove
- 2 Driver
- 3 Motor shaft
- 4 Hole

1. ➤ Open the lid.
2. ➤ Unscrew the rotor fastening nut using the supplied spanner.
3. ➤ Remove the rotor.

Installing the rotor

Personnel:

- Trained user

The lid is open.

1. ➤ Clean the motor shaft (3) and rotor hole (4).
2. ➤ Lightly grease the motor shaft (3), see ➔ Chapter 8.2 'Instructions for cleaning and disinfection' on page 25.
3. ➤ Place the rotor vertically on the motor shaft (3).
The driver (2) of the motor shaft must be in the groove (1) of the rotor.
4. ➤ Hand-tighten the rotor fastening nut using the supplied spanner.
5. ➤ Check that the rotor is firmly seated.

6.3 Loading

Filling centrifuge tubes

WARNING

Risk of injury from contaminated sample material.

Contaminated sample material escapes from the sample tube during centrifugation.

- Use centrifuge tubes with special screw caps for hazardous substances.
- For risk group 3 and 4 materials, use a biosafety system in addition to the sealable centrifuge tubes (see WHO's 'Laboratory Biosafety Manual').

NOTICE

Damage to the device due to highly corrosive substances.

Highly corrosive substances may impair the mechanical strength of rotors, buckets and accessories.

- Do not centrifuge highly corrosive substances.

Standard glass centrifuge tubes can be loaded up to RCF 4000 (DIN 58970 part 2).

Personnel:

- Trained user

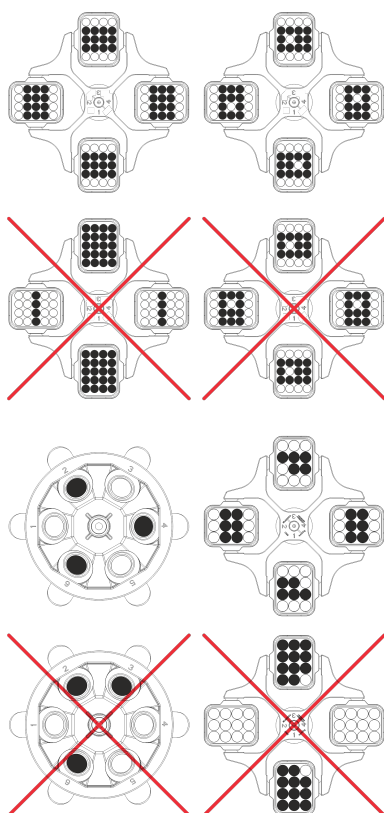
➤ Fill centrifuge tubes outside the centrifuge.

The maximum capacity of the centrifuge tubes specified by the manufacturer must not be exceeded.

With angle rotors, the centrifuge tubes must only be filled to the extent that no liquid can be ejected from the tubes during the centrifugation run.

It must be ensured that there is a uniform fill level in the tubes in order to keep the weight differences in the centrifuge tubes as low as possible.

Loading swing-out rotors

Personnel:

- Trained user

1. ➤ Check that the rotor is firmly seated.

2. ➤ The centrifuge tubes must be distributed symmetrically across all rotor locations.

The weight of the permissible filling capacity is indicated on each rotor. The weight must not be exceeded.

No liquid must be allowed to enter the buckets and the centrifuging chamber when loading the buckets and swinging them out during the centrifugation run.

For containers with rubber inserts, there must always be the same number of rubber inserts under the centrifuge tubes.

All rotor locations must be filled with the same buckets. Certain buckets are marked with the number of the rotor location. The buckets must only be inserted in the corresponding rotor location.

Buckets marked with a set number (for example S001/4) must only be used in the set.

6.4 Centrifugation

6.4.1 Centrifugation in continuous operation

Personnel:

- Trained user

1. ➤ Use the $[RPM \times 100]$ buttons to set the desired speed.

2. ➤ Use the $[t]$ buttons to set the time to zero.

➡ '0' is displayed.

3.  Press the *[START]* button.
 - The centrifugation run is started.

The 'Rotation' indicator light rotates while the rotor is turning.

Timing starts at 0. The first minute is counted up in seconds, then the time is displayed in minutes. A dot flashes next to the number if the time is displayed in minutes.

The rotor speed and the elapsed time are displayed during the centrifugation run.
4.  Press the *[STOP]* button to cancel the centrifugation run.

Ramp-down takes place with the set brake level.

An audible signal sounds when the rotor comes to a standstill.

6.4.2 Centrifugation with time preselection

Personnel:

- Trained user
1.  Use the *[RPM x 100]* buttons to set the desired speed.
 2.  Use the *[t]* buttons to set the desired time.

Adjustable from 1 to 99 minutes in 1 minute increments.
 3.  Press the *[START]* button.
 - The centrifugation run is started.

The 'Rotation' indicator light rotates while the rotor is turning.

The time is displayed in minutes. The last minute is counted down in seconds. A dot flashes next to the number if the time is displayed in minutes.

The rotor speed and the remaining time are displayed during the centrifugation run.
 4.  Ramp-down takes place with the selected brake level after the time has elapsed or if the centrifugation run is cancelled by pressing the *[STOP]* button.

An audible signal sounds when the rotor comes to a standstill.

6.4.3 Short-time centrifugation

Personnel:

- Trained user
1.  Use the *[RPM x 100]* buttons to set the desired speed.
 2.  Press and hold the *[PULSE]* button.
 - The centrifugation run is started.

The 'Rotation' indicator light rotates while the rotor is turning.

Timing starts at 0. The first minute is counted up in seconds, then the time is displayed in minutes. A dot flashes next to the number if the time is displayed in minutes.

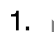
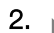
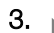
The rotor speed and the elapsed time are displayed during the centrifugation run.
 3.  Release the *[PULSE]* button to end the centrifugation run.
 - Ramp-down takes place with the set brake level. The brake level is displayed.

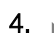
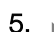
An audible signal sounds when the rotor comes to a standstill.

7 Software operation

7.1 Centrifugation parameters

7.1.1 Set the brake level

1.  Switch off the mains switch.
2.  Press and hold down the **▲** *[RPM x 100]* button and the *[PULSE]* button simultaneously.
3.  Switch on the mains switch and release the buttons.
 - Press the **▲** *[RPM x 100]* button repeatedly until the speed indicator shows the machine version (and/or 'C') and the time indicator shows the brake level set (and/or '0' or '1').

The machine version is set ex works and cannot be changed.
4.  Use the *[t]* buttons to set the desired brake level.
 - Level 1 = short ramp-down time.
 - Level 0 = long ramp-down time.
5.  Press the *[STOP]* button to save the settings.

7.1.2 Relative centrifugal force, RCF

The relative centrifugal force RCF is dependent on the speed and the centrifuging radius.

The relative centrifugal force RCF is stated as a multiple of the acceleration due to gravity (g).

The relative centrifugal force RCF is a dimensionless numerical value and is used to compare the separation and sedimentation performance.

$$RCF = \left(\frac{RPM}{1000} \right)^2 * r * 1,118$$

$$RPM = \sqrt{\frac{RCF}{r * 1,118} * 1000}$$

RCF = Relative Centrifugal Force

RPM = speed

r = centrifuging radius in mm = distance from the centre of the axis of rotation to the bottom of the centrifuge tube.

7.1.3 Centrifugation of substances or mixtures of substances with a density higher than 1.2 kg/dm³

The density of the substances or mixtures of substances must not exceed 1.2 kg/dm³ during centrifugation at maximum speed. The speed must be reduced for substances or substance mixtures with a higher density. The permissible speed can be calculated using the following formula:

$$\text{Reduced speed } (n_{red}) = \sqrt{\frac{1,2}{\text{Greater density [kg/dm}^3\text{]}} * \text{maximum speed [RPM]}}$$

For example: Maximum speed 4000 RPM, density 1.6 kg/dm³

$$n_{red} = \sqrt{\frac{1,2(\text{kg/dm}^3)}{1,6(\text{kg/dm}^3)} * 4000 \text{ RPM}} = 3464 \text{ RPM}$$

If, in exceptional cases, the maximum load indicated on the bucket is exceeded, the speed must also be reduced. The permissible speed can be calculated using the following formula:

$$\text{Reduced speed } (n_{red}) = \sqrt{\frac{\text{maximum load [g]}}{\text{actual load [g]}}} * \text{maximum speed [RPM]}$$

For example: Maximum speed 4000 RPM, maximum load 300 g, actual load 350 g

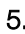
$$n_{red} = \sqrt{\frac{300 \text{ g}}{350 \text{ g}}} * 4000 \text{ RPM} = 3703 \text{ RPM}$$

Please contact the manufacturer if you are not sure.

7.1.4 Centrifuging radius

The centrifuging radius must be entered in centimetres.

It is possible to enter values between 5 cm and 16 cm.

1.  Switch off the mains switch.
2.  Press and hold down the  *[RPM x 100]* button and the *[PULSE]* button simultaneously.
3.  Switch on the mains switch and release the buttons.
4.  Press the  *[RPM x 100]* button repeatedly until the centrifuging radius and 'rd' are displayed.
5.  Use the *[t]* buttons to set the desired centrifuging radius.
6.  Press the *[STOP]* button to save the settings.

7.2 Machine Menu

7.2.1 Audible signal

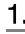
7.2.1.1 General

The audible signal sounds:

- after a problem occurs in the 2 s interval.
- after completion of the centrifugation run and rotor standstill in the 30 s interval.

Opening the lid or pressing any button stops the audible signal.

7.2.1.2 Setting an audible signal

1.  Switch off the mains switch.
2.  Press and hold down the  *[RPM x 100]* button and the *[PULSE]* button simultaneously.
3.  Switch on the mains switch and release the buttons.
4.  Press the  *[RPM x 100]* button repeatedly until 'BEL 1' or 'BEL 0' is displayed.
5.  Set '0' or '1' using the *[t]* buttons beneath the time indicator.
0 = audible signal disabled.
1 = audible signal enabled.
6.  Press the *[STOP]* button to save the settings.

8 Cleaning and care

8.1 Overview table

| Chap. | Task to execute | if required | daily | weekly | Annually | Page |
|------------|--|-------------|-------|--------|----------|------|
| 8 | Cleaning and care | | | | | 25 |
| 8.3 | Cleaning | | | | | 26 |
| 8.3 | Cleaning the device | | X | | | 26 |
| 8.3 | Cleaning accessories | | | X | | 26 |
| 8.4 | Disinfection | | | | | 26 |
| 8.4 | Disinfecting the device | X | | | | 27 |
| 8.4 | Disinfecting the accessories | X | | | | 27 |
| 8.5 | Maintenance | | | | | 27 |
| 8.5 | Greasing the rubber seal of the centrifuging chamber | | | X | | 27 |
| 8.5 | Trunnion greasing | | | X | | 27 |
| 8.5 | Checking the accessories | | | X | | 27 |
| 8.5 | Centrifuging chamber damage inspection | | | | X | 27 |
| 8.5 | Greasing the motor shaft | | | | X | 27 |
| 8.5 | Accessories with a limited service life | X | | | | 27 |
| 8.5 | Replacing centrifuge tubes | X | | | | 28 |

8.2 Instructions for cleaning and disinfection

DANGER

Risk of contamination for the user due to inadequate cleaning or failure to observe the cleaning instructions.

- Observe cleaning instructions.
- Wear personal protective equipment when cleaning the device.
- Observe laboratory regulations (e.g. TRBAs, the German Protection against Infection Act, hygiene plan) for handling biological agents.

- The device and its accessories must not be cleaned in dishwashers.
- Only perform hand cleaning and liquid disinfection.

- The water temperature must not exceed 25 °C.
- To prevent any corrosion due to use of detergents or disinfectants, it is essential to follow the special application instructions provided by the manufacturers of the detergent or disinfectant.

Disinfectant:

- Surface disinfectant (not disinfectant for hands or instruments)
- Ethanol as the sole active substance.
Do not use an ethanol-propanol mixture to disinfect the viewing window in the lid of the device.
- Concentration not less than 30%
- pH: 6 – 8
- Non-corrosive

8.3 Cleaning

Cleaning the device

1. ➤ Open the lid.
2. ➤ Switch off the device and disconnect it from the power supply.
3. ➤ Remove accessories.
4. ➤ Clean the centrifuge housing and the centrifuging chamber with soap or a mild detergent and a damp cloth.
5. ➤ Remove any detergent residues with a damp cloth after using detergents.
6. ➤ The surfaces must be dried immediately after cleaning.
7. ➤ Dry the centrifuging chamber with an absorbent cloth if condensation forms.

Cleaning accessories

1. ➤ Clean the accessories using the detergent and a damp cloth.
2. ➤ Remove any detergent residues with a damp cloth after using detergents.
3. ➤ Dry the accessories immediately after cleaning using a lint-free cloth and oil-free compressed air. Dry all cavities completely using oil-free compressed air.

8.4 Disinfection

Disinfection must always be preceded by cleaning the components concerned.

See ➔ Chapter 8.3 'Cleaning' on page 26

Disinfectant concentration and application time according to the manufacturer's instructions.

Disinfecting the device**CAUTION****Risk of injury due to ingress of water or other liquids.**

- Protect the device against external liquids.
- Do not disinfect the device using spray.

1. Open the lid.
2. Switch off the device and disconnect it from the power supply.
3. Remove accessories.
4. Clean the housing and centrifuging chamber using disinfectant.
5. Remove any disinfectant residues with a damp cloth after using disinfectants.
6. The surfaces must be dried immediately after cleaning.

Disinfecting the accessories

1. Disinfect the accessories using the disinfectant.
2. Wet all cavities with bubble-free disinfectant.
3. Remove the disinfectant residues or leave them to dry after using disinfectants.

Autoclaving

The rotor and accessories must not be autoclaved.

8.5 Maintenance**Greasing the rubber seal of the centrifuging chamber**

- Rub the sealing ring lightly with a rubber care product.

Trunnion greasing

1. Remove accessories.
2. Clean the trunnions.
3. Remove any detergent residues with a damp cloth after using detergents.
4. Grease the trunnions and suspension with Hettich Tubenfett 4051.
5. Excess grease in the centrifuging chamber must be removed.

Checking the accessories

1. The accessories shall be checked for wear and corrosion damage.
2. Check that the rotor is firmly seated.

Centrifuging chamber damage inspection

- Check the centrifuging chamber for damage.

Greasing the motor shaft

1. Remove accessories.
2. Clean the motor shaft.
3. Remove any detergent residues with a damp cloth after using detergents.
4. Grease the motor shaft with Hettich Tubenfett 4051.
5. Excess grease in the centrifuging chamber must be removed.

Accessories with a limited service life

The use of certain accessories is time-limited. For safety reasons, the accessories must no longer be used when either the maximum number of permissible run cycles marked on them or the expiry date marked on them has been reached.

- The maximum permissible number of run cycles or the expiry date can be seen on the accessories.
- The centrifuge is equipped with a cycle counter.

Replacing centrifuge tubes

CAUTION

Risk of injury from broken glass.

Broken glass may cause glass splinters and contaminated liquids to be found inside the centrifuge.

- Wear cut-resistant gloves.
- Wear protective goggles and a face mask.

Broken parts of the tube, glass splinters and spilled centrifuge material must be removed completely in the event of leakage or if a centrifuge tube breaks. Glass splinters that are not removed will cause further glass breakage.

The rubber inserts and the plastic sleeves of the rotors must be replaced after a glass breakage.

Disinfection must be carried out if the material is infectious.

9 Troubleshooting

9.1 Fault description

Customer service must be notified if the fault cannot be rectified based on the fault table. State the centrifuge type and serial number. Both numbers can be seen on the type plate of the centrifuge.

| Fault description | Cause | Remedy |
|-------------------|--|---|
| no display | No power. Mains input fuses defective. | <ul style="list-style-type: none"> ■ Check the supply voltage. ■ Check the mains input fuse. ■ Set the mains switch to <i>/I/</i>. |
| -1- | Tacho error. Speed pulse failure. | <ul style="list-style-type: none"> ■ The device must not be switched off while the 'Rotation' indicator is lit up and rotating. <p>Wait until the 'Lid closed' symbol is displayed (after approx. 120 seconds) and then perform a MAINS RESET.</p> |
| -2- | Loss of mains power during the centrifugation run. The centrifugation run was not completed. | <ul style="list-style-type: none"> ■ Open the lid and press the <i>[START]</i> button. ■ If required: Repeat the centrifugation run. |
| -3- | Imbalance. The rotor is unevenly loaded. | <ul style="list-style-type: none"> ■ Open the lid. ■ Check the loading of the rotor. ■ Repeat the centrifugation run. |
| -4- | Communication. Error in the control section or power section. | <ul style="list-style-type: none"> ■ Perform a MAINS RESET. |
| -5- | Overload. Motor or motor control defective. | <ul style="list-style-type: none"> ■ Perform a MAINS RESET. |

| Fault description | Cause | Remedy |
|--|---|---|
| -6- | Overvoltage. Mains voltage outside tolerances. | <ul style="list-style-type: none"> ■ Perform a MAINS RESET. ■ Check the mains voltage. |
| -7- | Overspeed. Error in the power section. | <ul style="list-style-type: none"> ■ Perform a MAINS RESET. |
| -8- | Undervoltage. Mains voltage outside tolerances. | <ul style="list-style-type: none"> ■ Perform a MAINS RESET. ■ Check the mains voltage. |
| -9- | Overtemperature. Overtemperature switch in the motor has tripped. | <ul style="list-style-type: none"> ■ Open the lid using the emergency release. ■ Let the motor cool down. |
| Version Error | Wrong machine version set. The control section skips to the Settings menu. | <ul style="list-style-type: none"> ■ Use the <i>[t]</i> buttons to set the letter C. ■ Press the <i>[STOP]</i> button to save the settings. ■ Perform a MAINS RESET. |
| no speed indicator. Machine version set in the time indicator. | Version Error. Wrong machine version set. The control section skips to the Settings menu. | <ul style="list-style-type: none"> ■ Use the <i>[t]</i> buttons to set the letter C. ■ Press the <i>[STOP]</i> button to save the settings. ■ Perform a MAINS RESET. |
| -b- | Underspeed. | <ul style="list-style-type: none"> ■ Perform a MAINS RESET. |
| -c- | Controller watchdog. Error in the power section. | <ul style="list-style-type: none"> ■ Perform a MAINS RESET. |
| -d- | Lid lock error. | <ul style="list-style-type: none"> ■ Perform a MAINS RESET. |
| -E- | Short circuit in the control section / power section. | <ul style="list-style-type: none"> ■ Perform a MAINS RESET. |
| -F- | Wrong machine version. | <ul style="list-style-type: none"> ■ Notify customer service. |
| 888888 All indicator segments light up. | - | <ul style="list-style-type: none"> ■ Notify customer service. |

9.2 Performing a MAINS RESET

1. ➤ Set the mains switch to *[0]*.
2. ➤ Wait 10 seconds.
3. ➤ Set the mains switch to *[I]*.

9.3 Emergency release

The lid cannot be unlocked by the motor in the event of a power failure. Emergency unlocking by hand must be performed.

WARNING

Risk of electric shock due to maintenance and servicing work on live device.

- Disconnect the device from the mains before carrying out repairs and maintenance.

WARNING

Danger of cutting and crushing due to moving rotor.

- Do not open the lid until the rotor has stopped.

The plastic unlocking pin supplied may only be used for emergency device release. The release pin must be stored in such a way that it is protected against unauthorised access.

Personnel:

- Trained user

1. Look through the window in the lid to ensure that the rotor is stationary.
2. Insert the supplied plastic release pin horizontally into the hole (1). Push it in until the lid can be opened when the pin is pressed down.
3. Open the lid.

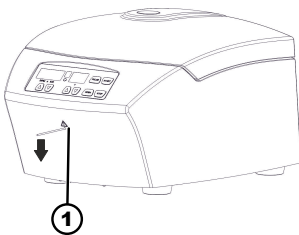
Fig. 16: Emergency release

1 Hole

9.4 Replacing the mains input fuse

WARNING

Risk of electric shock due to maintenance and servicing work on live device.

- Disconnect the device from the mains before carrying out repairs and maintenance.

Personnel:

- Trained user

The mains fuses are located next to the mains switch.

The mains switch is in switch position [O]

1. Disconnect the mains cable from the device plug.
2. Press the snap locks (2) against the fuse holder (1) and pull them out.
3. Replace the defective mains input fuses.
Only use fuses with the nominal value specified for the type: see the table below.
4. Push in the fuse holder (1) until the snap lock engages.
5. Reconnect the device to the mains.

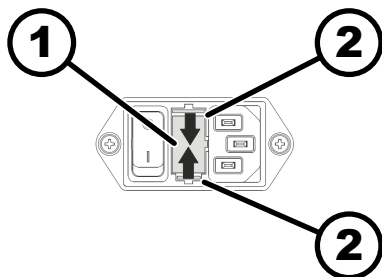
Fig. 17: Mains input fuse

1 Fuse holder
2 Snap lock

| Model | Type | Fuse | Order no. |
|---------|------|----------------|-----------|
| EBA 270 | 2300 | T 1.6 AH/250 V | E891 |

| Model | Type | Fuse | Order no. |
|---------|---------|--------------------|-----------|
| EBA 270 | 2300-01 | T 3.15 AH/250 V | E997 |

10 Disposal

10.1 General instructions

The device can be disposed of via the manufacturer.

A Return Material Authorisation (RMA) form must always be requested for a return.

If necessary, contact the Technical Service Department of the manufacturer:

- **Andreas Hettich GmbH & Co. KG**
- Föhrenstraße 12
- 78532 Tuttlingen, Germany
- Phone: +49 7461 705 1400
- E-Mail: service@hettichlab.com

! WARNING

Risk of pollution and contamination for people and the environment.

When disposing of the centrifuge, people and the environment may be polluted or contaminated by incorrect or improper disposal.

- Removal and disposal may be carried out only by a trained and authorized service personnel.

The device is intended for the commercial sector ("Business to Business" - B2B).

According to Directive 2012/19/EU, the devices may no longer be disposed of with household waste.

The appliances are assigned to the following groups according to the Stiftung Elektro-Altgeräte Register (EAR (German foundation under civil law)):

- Group 5 (small appliances)

The crossed-out wheelee bin symbol indicates that the device must not be disposed of with household waste. Regulations governing disposal of such devices may differ in individual countries. If necessary, contact the supplier.

Fig. 18: Household waste ban

11 Index

| | | | |
|--|----|--|--------|
| A | | M | |
| Accessories. | 14 | MAINS RESET. | 29 |
| biosafety systems. | 26 | Maintenance. | 27 |
| disinfection. | 27 | Intervals. | 25 |
| the rotor. | 27 | N | |
| with limited service life. | 27 | Non-intended use. | 6 |
| Autoclaving. | 27 | O | |
| C | | Operator's responsibility. | 7 |
| Care | | Original spare parts. | 14 |
| Intervals. | 25 | P | |
| Centrifugation | | Personal protective equipment. | 6 |
| in continuous operation. | 21 | Personnel qualifications. | 6 |
| with higher substance density. | 23 | Personnel training. | 7 |
| with time preselection. | 22 | Protective equipment. | 6 |
| Centrifuging chamber | | R | |
| inspection. | 27 | Rating plate. | 10 |
| Cleaning. | 26 | Relative centrifugal force | |
| Cleaning and disinfection | | RCF. | 23 |
| Instructions. | 25 | Removing | |
| Connecting the centrifuge. | 18 | the rotor. | 19, 21 |
| Continuous operation. | 21 | Replacing | |
| D | | centrifuge tubes. | 28 |
| Device | | Return. | 14 |
| biosafety systems. | 26 | S | |
| disinfection. | 27 | Safety instructions. | 7 |
| Disinfection. | 26 | Scope of delivery. | 14 |
| Disposal. | 31 | Setting up the centrifuge. | 17 |
| E | | Short-term centrifugation. | 22 |
| Enabling/disabling. | 24 | Spare parts. | 14 |
| Error messages. | 28 | Storage conditions. | 15 |
| F | | Switching off. | 19 |
| Fastening | | Switching the unit on. | 19 |
| removal. | 17 | Symbols. | 5 |
| the transport lock. | 15 | T | |
| Filling. | 20 | Transport condition. | 15 |
| Foreseeable misuse. | 6 | Troubleshooting. | 28 |
| G | | Trunnion | |
| General safety instructions. | 7 | greasing. | 27 |
| Greasing | | U | |
| the rubber seal. | 27 | Unpacking. | 16 |
| I | | | |
| Intended use. | 5 | | |
| L | | | |
| Labels | | | |
| on the device. | 12 | | |
| on the packaging. | 11 | | |
| Lid | | | |
| the lid. | 19 | | |
| Loading. | 20 | | |

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