MIKRO 2.0 / 2.0 R

# Inhalt des Dokuments / content of the document

Gebrauchsanweisung (DE) Operating instructions (EN) Mode d'emploi (FR) Istruzioni per l'uso (IT) Rotoren und Zubehör / Rotors and accessories

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MIKRO 2.0 / 2.0 R

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# Operating instructions MIKRO 2.0 / 2.0 R

Translation of the original operating instructions

©2025 - All rights reserved Andreas Hettich GmbH Föhrenstrasse 12 D-78532 Tuttlingen, Germany Telephone: +49 (0)7461 705-0 Fax: +49 (0)7461 705-1125 Email: info@hettichlab.com, service@hettichlab.com Internet: www.hettichlab.com

Table of contents

1	А	bout this document	5
	1.1	Use of this document	5
	1.2	Gender reference	5
	1.3	Symbols and labels in this document	5
	1.4	Explanation of terms	6
2	S	afety	6
	2.1	Intended use	6
	2.2	Foreseeable misuse	6
	2.3	Use of accessories	7
	2.4	Service life	7
	2.5	Personnel qualifications	7
	2.	5.1 Authorisation levels	7
	2.6	Personal protective equipment.	8
	2.7	Operator obligations	8
	2.8	Safety instructions	9
3	D		11
	3.1	Technical data	11
	3.2	Rating plate	12
	3.3	Certifications and logos.	13
	3.4	European registration	14
	3.5	Important labels on the packaging	14
	3.6	Important labels on the device	15
	3.7	Returns	15
	3.8	Operating and indicator elements.	16
	3.	8.1 Control	16
	3.	8.2 Controls	16
	3.9		17
4	TI 		17
	4.1		17
	4.2		18
_	4.3		18
5			19
	5.1		19
	5.2		19
	5.3	Setting up and connecting the centrifuge	19
	5.4	Switching on for the lifst time	21
e	0.0		22
Ö		Opening and electing the lid	23
	0.1		23
	0.2		24 05
	0.3 6 1		20 20
	0.4		20

6	.5 Cer	ntrifugation parameters	27
	6.5.1	Relative centrifugal force, RCF	27
	6.5.2	Centrifugation of substances or mixtures of substances	
		with a density higher than 1.2 kg/dm <sup>3</sup>	28
	6.5.3	Ramp-up and brake levels	29
6	.6 Use	er interface	29
	6.6.1	Dealing with messages	29
	6.6.2	Basic displays	30
	6.6.3	Keypad	32
6	.7 Mer	nus	33
	6.7.1	Main menu	33
	6.7.2	Quick Adjust menu	34
	6.7.3	Program menu	34
	6.7.4	Settings menu	35
6	.8 Cer	ntrifugation	35
	6.8.1	Continuous operation	35
	6.8.2	Centrifugation with time preselection	36
	6.8.3	Short-term centrifugation	37
6	.9 Qui	ck stop function	38
6	.10 Soff	tware operation	39
	6.10.1	Set Quick Adjust	39
	6.10.2	Program menu for authorisation levels	41
	6.10.3	Device information.	45
6	.11 Use	er settings	47
6	.12 Ope	en source licences	58
7	Mainte	enance, cleaning and care	58
7	.1 Clea	aning	59
7	.2 Disi	nfection	59
7	.3 Mai	ntenance	61
8	Faults		63
8	.1 Fau	It description	63
8	.2 War	rning descriptions	67
8	.3 Per	forming a MAINS RESET	71
8	.4 Eme	ergency release	71
8	.5 Rep	placing the mains input fuse	72
9	Disass	sembly	72
10	Spare	parts	72
11	Stora	ge	73
12	Dispo	sal	73
1	2.1 Ger	neral instructions	73
13	Gloss	ary	75
14	Index.		76

## 1 About this document

## 1.1 Use of this document

- Read this document carefully and in full before commissioning the device for the first time.
  - Observe other enclosed instruction sheets where necessary.
- This document is part of the device and must be kept within easy reach.
- This document must be included if the device is passed on to a third party.
- The current version of the document in the available languages can be found on the manufacturer's website: *https://www.hettichlab.com/de/ download-center/*

## 1.2 Gender reference

The employed masculine or feminine language form is to facilitate reading. In the spirit of equal treatment, corresponding terms apply in principle to all genders and do not imply any valuation.

### 1.3 Symbols and labels in this document

#### General symbols

The following markers are used in this document to highlight instructions, results, listings, references and other elements:

Marker	Explanation
1 2 3	Step-by-step instructions
⇒	Results of action steps
►	References to sections of the document and other applicable documents
•	Listings without a fixed order
[Buttons]	Controls (for example: buttons, switches)
'Indicator'	Indicator elements (for example: signal lights, screen elements)

#### Warnings

Hazard levels	Meaning
DANGER	Indicates a hazard which, if not avoided, will result in death or serious injury.
WARNING	Indicates a hazard which, if not avoided, could result in death or serious injury.
CAUTION	Indicates a hazard which, if not avoided, could result in minor or moderate injuries.

Hazard levels	Meaning
NOTE	Indicates a hazard which, if not avoided, could result in damage to property.

#### Worth knowing

Recommenda- tions	Meaning
Tip/Info	Indicates information for efficient operation and know-how for effective use of the device.

## 1.4 Explanation of terms

The term "centrifuge" employed in these operating instructions refers to the device referred to on the cover sheet.

## 2 Safety

2.1 Intended use

The centrifuge **MIKRO 2.0 General Purpose**, **MIKRO 2.0 R General Purpose** is a general laboratory device that is placed on the market in accordance with Machinery Directive 2006/42/EC (a changeover to Machinery Directive 2023/1230 will occur in July 2026).

The device is used to separate aqueous solutions and suspensions up to a density of 1.2 kg/dm<sup>3</sup>. The user can set the variable physical parameters within the limits specified by the device.

The centrifuge may only be used by qualified personnel in closed laboratories.

**Disclaimer** Intended use also includes observing all information in the operating instructions and performance of all the required inspection and maintenance work. Any other use or use beyond this scope is considered improper. Andreas Hettich GmbH shall not be liable for any damage arising from such noncompliant use.

If the centrifuge is installed in another device or integrated into a system, the manufacturer of the system as a whole is responsible for its safety.

## 2.2 Foreseeable misuse

Use of the centrifuge outside the defined conditions is considered improper use.

Among other things, the following are considered improper use:

- Use of the centrifuge in explosive environments.
- Use of the centrifuge in radioactive environments.
- Use of the centrifuge in biologically contaminated environments.
- Use of the centrifuge in chemically contaminated environments.
- Use of the centrifuge in corrosive environments.
- Centrifugation of combustible materials.
- Centrifugation of explosive materials.
- Centrifugation of radioactive materials.
- Centrifugation of chemically highly reactive materials.
- Centrifugation of corrosive materials.

2.3	Use of accessories	
		Only use accessories approved by the manufacturer for the intended use of the centrifuge.
2.4	Service life	
		The planned service life of the device is 7 years.
		Service life of the accessory in terms of the number of cycles and/or service life as indicated on the accessory.
2.5	Personnel qualificati	ons
		All control and safety devices must only be operated by trained persons.
	1	
Qualifie	d mechanics	For the purposes of this documentation, qualified mechanics are persons who are familiar with the design, mechanical installation, commissioning, troubleshooting and maintenance of the product and who have the fol- lowing qualifications:
		<ul> <li>Qualification / training in the field of mechanics in accordance with national regulations.</li> </ul>
Qualifie	d electricians	For the purposes of this documentation, qualified electricians are persons who are familiar with the design, electrical installation, commissioning, trou- bleshooting and maintenance of the product and who have the following qualifications:
		<ul> <li>Qualification / training in electrical engineering in accordance with national regulations.</li> </ul>
Qualifie cians	d refrigeration techni-	For the purposes of this documentation, qualified refrigeration technicians are persons who are familiar with the design and installation of refrigeration and chilling equipment, and the commissioning, troubleshooting and main- tenance of the product and who have the following qualifications:
		<ul> <li>Qualification / training in the field of refrigeration and cooling technology in accordance with national regulations.</li> </ul>
		<ul> <li>An appropriate certificate of competence (e.g. in accordance with ISO 22712) for working with A3 refrigerants must be available.</li> </ul>
Trained	person	For the purposes of this documentation, these are persons who have received instruction enabling them to carry out work in the areas of trans- portation, storage and operation.
2.5.1	Authorisation levels	
Authori	sation level 1	"Restricted rights"
		A user who uses the system to complete a task. Centrifugations can be started and stopped. Changing programs or viewing and changing system settings is not permitted.
		The user can only use some functions after a user with authorisation level 3 has entered a PIN.
		Example: loader / assistant / student

Authorisation level 2 "Standard rights"

Authorisation level 1 rights are included.

A user with advanced knowledge. The user is allowed to change the value settings in Quick Adjust. This has a corresponding effect on short-term centrifugation. Event messages/logs can be viewed. Readout and analysis of system information is possible to a limited extent.

Some functions can only be used by authorisation level 2 users after an authorisation level 3 user has entered a PIN.

Authorisation level 3 Authorisation level 2 rights are included.

A user who is allowed to make changes to the system and programs. Readout and analysis of system information is possible.

The user must enter a PIN on initial start-up.

Example: laboratory manager / product manager

## 2.6 Personal protective equipment

Observe national and local regulations on safety and accident prevention. Select and provide protective clothing such as protective footwear, safety gloves and protective goggles in alignment with the respective activities and the expected risks.

### 2.7 Operator obligations

Ensure that all the work listed below is carried out only by qualified personnel:

- Transportation, unpacking, lifting
- Setup
- Operation
- Maintenance
- Disposal

The operator must ensure that persons working on the device observe the rules and regulations as well as the following instructions:

- National and local regulations pertaining to safety, accident prevention and environmental protection.
- Do not install or commission any damaged devices.
- The necessary protective equipment must be provided.
- Operate the device only when it is in perfect working order.
- Periodically check that personnel are working safely and in compliance with the safety information in the operating instructions.
- Check protective equipment regularly to ensure it is serviceable.
- Do not remove attached safety and warning notices and ensure they remain legible.
- Persons under the influence of alcohol, drugs or medicines that affect their ability to react must not operate or maintain the device.
- Keep the operating instructions available in the place where the device is used.
- Pass the operating instructions on to the buyer when the device is sold.

#### Reporting of events and incidents

In the event of incidents or notifiable events involving the device or its accessories, these must be reported to the manufacturer and, where applicable, to the competent authority where the user and/or the patient is registered. Manufacturer:

Safety

Andreas Hettich GmbH

Föhrenstrasse 12 78532 Tuttlingen, Germany Phone: +49 7461 705 0 E-mail: info@hettichlab.com

#### DANGER

#### Risk of contamination

Inadequate cleaning or failure to observe the cleaning instructions can lead to contamination risks.

- Observe national and local regulations on safety and accident prevention.
- Observe cleaning instructions.
- Wear personal protective equipment when cleaning the device.
- Comply with laboratory regulations (e.g. TRBAs, the German Protection against Infection Act, hygiene plan) for handling biological agents.

## DANGER

#### Risk of fire and explosion

Risk of accidents, injuries or damage to property due to fire or explosion.

- Observe regulations and directives for handling chemicals and hazardous substances.
- Do not use corrosive chemicals.
- Do not use dangerous chemicals.
- Do not use corrosive extraction agents.
- Do not use strong acids.

# WARNING

## Risk of injury

Inadequate or late maintenance can result in injuries.

- Observe maintenance intervals.
- Check the device for visible damage or defects.
  - If any visible damage or defects are present, immediately remove the device from service and inform a service technician.

## 🕂 WARNING

#### Risk of electric shock

Liquids that penetrate the device may cause electric shocks.

- The device must be protected from external contact with liquids.
- Do not pour any liquids inside the device.
- The original transport packaging must be used when transporting the device.

## WARNING

# Contamination with hazardous substances and substance mixtures

Observe the following actions for substances and substance mixtures that are toxic, radioactive and/or contaminated with pathogenic microorganisms:

- As a rule, use only centrifuge tubes with special screw caps for hazardous substances.
- Use sealable centrifuge tubes with a biosafety system for materials of risk groups 3 and 4.
- If no biosafety system is used, the device is not microbiologically tight in the sense of standard EN / IEC 61010-2-020.
- Contact the manufacturer if necessary.

## Risk of injury and device damage due to the rotor

A loose rotor can cause injuries and device damage.

- Check that the rotor is firmly seated.
- Follow maintenance intervals.

## CAUTION

#### Risk of injury

Long hair and clothing can get caught in the rotor during manual movement.

- Tie long hair back.
- Do not allow garments to hang in the centrifuging chamber.

#### NOTICE

#### Damage

Incorrect voltage or frequency.

 Only operate the device according to the specifications on the rating plate.

Compliance with the instructions for use.

NOTICE

## Damage

Aborting the program early can cause damage to the device and samples.

- Do not switch off, perform an emergency release or pull out the mains plug.

## 3 Device overview

# 3.1 Technical data

Manufacturer	Andreas Hettich GmbH D-78532 Tuttlingen		
Model	MIKRO 2.0	MIKRO 2.0 R	
Туре	2440	2460	
Mains voltage (±10%)	200-240 V 1~	200-240 V 1~	
Mains frequency	50-60 Hz	50-60 Hz	
power consumption	270 VA	420 VA	
Refrigerant	-	R290	
Refrigerant content	-	35 g	
Global warming potential (GWP)	-	0,02	
max. capacity	12 x 5 ml		
max. permissible density	1.2 kg/dm <sup>3</sup>		
max. speed (RPM)	16100		
max. acceleration (RCF)	25000		
max. kinetic energy	6400 Nm		
Obligation to perform checks (DGUV Rules 100-500) (valid only in Germany)	No		
Ambient conditions (EN / IEC 61010-1):			
Installation site	indoors only		
Altitude	up to 2000 m above sea level		
Ambient temperature	2 °C to 40 °C	5 °C to 35 °C	
Humidity	maximum relative humidity 80% for temperatures up to 31 °C, decreasing linearly to 50% relative humidity at 40 °C.		

Overvoltage category (IEC 60364-4-443)	II	
Pollution level	2	
Device protection class	I not suitable for use in potentially explosive atmospheres.	
EMC:		
Emitted EM interference, EM interference immunity	EN / IEC 61326-1 Class B FCC Class B	EN / IEC 61326-1 Class B FCC Class B
Noise level (rotor-dependent)	≤58 dB(A)	≤61 dB(A)
Dimensions:		
Width	240 mm	240 mm
Depth	389 mm	538 mm
Altitude	225 mm	250 mm
Weight	approx. 18 kg	approx. 28 kg

# 3.2 Rating plate

Fig. 1: Rating plate

1	REF Item number	12	Maximum permissible density
2	Serial number	13	Manufacturer's address
3	Revision	14	[]] Follow the user manual
4	Equipment number	15	Maximum permissible pressure (PS) in the refrigeration system (only for devices with cooling)

5	UDI code	16	Refrigerant content (only for devices with cooling)
6	Content of the UDI code: (01) Global Trade Item Number (GTIN) (11) Date of manufacture (21) Serial number	17	Refrigerant type (only for devices with cooling)
7	EAC mark, CE mark	18	Maximum speed in revs per minute
8	Country of manufacture	19	Current draw (A) or power consumption (VA)
9	<sup>™</sup> Date of manufacture	20	Device designation
10	Mains voltage and mains frequency	21	Manufacturer's logo
11	Maximum kinetic energy		

# 3.3 Certifications and logos

## Certifications

G TUV SUD ISO 9001	O 9001 tified ality Management System w.tuvsud.com/ms-cert	ISO 9001 Quality management system in accordance with ISO 9001
ISO ISO 14001	O 14001 tified Environmental hagement System w.tuvsud.com/ms-cert	ISO 14001 Environmental management in accordance with ISO 14001
Certified EN ISC Quality System Quality	) 13485 v management in accorda	nce with ISO 13485

#### Logos

**Omdc** 

## 3.4 European registration

**Device conformity** Device conformity according to EU directives.

TOP

## 3.5 Important labels on the packaging

This is the correct upright position of the shipping container for transport and/or storage.

FRAGILE GOODS The contents of the shipping container are fragile, so it must be handled with care.

### PROTECT FROM MOISTURE

Protect the transport packaging from moisture and keep it in a dry environment.

#### TEMPERATURE LIMITATION

The shipping container must be stored, transported and handled within the indicated temperature range (-20  $^\circ C$  to +60  $^\circ C$ ).

#### HUMIDITY LIMITATION

The shipping container must be stored, transported and handled within the indicated air humidity range (10 % to 80 %, non-condensing).

#### STACK LIMITATION BASED ON QUANTITY

Maximum number of identical packages that may be stacked on the lowest package, "n" standing for the number of packages allowed. The lowest package is not included in "n".

Warning of risk of fire and explosion due to flammable refrigerant

## 3.6 Important labels on the device

(	$\bigcirc$	)
_		

The signs on the device must not be removed or covered, or have anything pasted over them.

Attention, general danger area.

Ensure you read the instructions for commissioning and operation and observe the safety instructions before using the device.

Biohazard warning.

Warning of fire hazard.

This warning indicates the risk of fire or explosion caused by refrigerants.

Direction of rotation of the rotor.

The orientation of the arrow indicates the rotor's direction of rotation.

Direction of rotation of the emergency release.

Symbol for the separate collection of electrical and electronic equipment, in accordance with Directive 2012/19/EU (WEEE).

Use in European Union countries, Norway and Switzerland.

## 3.7 Returns

An original Return Material Authorisation (RMA) form from the manufacturer must always be requested for a return. Secure and reliable acceptance and booking in of the goods with the manufacturer is not possible without an original RMA form from the manufacturer. The Return Material Authorisation (RMA) form contains a Declaration of No Objection (UBE), which must be completed in full and enclosed with the return.

If a leak is suspected or confirmed by the service department, the refrigeration circuit must be evacuated prior to return. The same applies in the event of a crash because it is not certain that all the refrigerant will escape.

If the device and/or accessories are returned to the manufacturer, the complete return shipment must be cleaned and decontaminated by the sender. If returns are not cleaned and/or decontaminated or are insufficiently cleaned and/or decontaminated, this will be performed by the manufacturer and charged to the sender.

The original transport locks must be attached for return shipment, see → Chapter 4 'Transport and storage' on page 17. The device must be shipped in its original packaging.

3.8 Operating and indicator elements

### 3.8.1 Control

Fig. 3: Control (device with cooling)

## 3.8.2 Controls

Number	Button and display ele- ments	Description
1	Screen	<ul> <li>Displays, information and menus</li> </ul>
2	Rotary and push button	<ul><li>Rotate: set the parameters and change to the next menu item.</li><li>Press: jump to the next parameter and select a menu item.</li></ul>

Number	Button and display ele- ments	Description
3	STOP, OPEN	<ul> <li>Press once: centrifuge brakes with set braking ramp.</li> <li>double or long press (quick stop): centrifuge brakes with maximum braking ramp.</li> <li>Press when the centrifuge is at a standstill: open the lid.</li> </ul>
4	START	<ul> <li>Start centrifugation run.</li> <li>A centrifugation run can be started from Quick Adjust or from the program.</li> </ul>
5	SHORT-TERM CENTRIFU- GATION	<ul> <li>The centrifugation run takes place as long as the button is being pressed.</li> </ul>
6	BACK	<ul><li>Back to the menus.</li><li>Save changed values.</li></ul>
7	INFO	<ul> <li>Operational data and rotor data retrieved.</li> </ul>
8	CONTEXT BUTTON, TOGGLE BUTTON	<ul> <li>You can access the context menu by pressing the button as soon as '' is displayed in a menu.</li> <li>To select the second level of favourite programs, hold down the button and simultaneously press the desired program shortcut button.</li> </ul>
9	SHORTCUT BUTTONS	<ul> <li>Shortcut buttons for favourite programs P1-P8.</li> <li>Shortcut buttons for favourite programs P1-P7 for MIKRO 2.0 R only.</li> <li>Temperature control program for MIKRO 2.0 R only.</li> </ul>
10	QUICK ADJUST, PRO- GRAM	<ul> <li>Quick Adjust: Quick setting of parameters. The values can subsequently be transferred to a program.</li> <li>Program: Setting parameters that are saved as a program. The program must be saved with a program name. Programs can be saved as favourites.</li> </ul>

## 3.9 Scope of delivery

- 1 offset screwdriver
- 1 power cable
- 1 user manual
- For products distributed in Switzerland:
   1 Import Switzerland sticker

# 4 Transport and storage

4.1 Transport and storage conditions

Transport conditions

#### NOTICE

The centrifuge's refrigeration circuit contains flammable refrigerant.

- Comply with possible transport restrictions (e.g. according to ADR or IATA DGR).

	NOTICE
	Damage
	<ul> <li>The device may be damaged if it is not secured during transport.</li> </ul>
	<ul> <li>Secure the transport locks before transport.</li> </ul>
	<ul> <li>Observe the transport instructions.</li> </ul>
	NOTICE
	<ul> <li>Danger due to condensation in the event of temperature differences</li> </ul>
	Moisture may damage electrical components.
	<ul> <li>Ensure that all surfaces are dry prior to commissioning or maintenance.</li> </ul>
	<ul> <li>If the temperature changes, wait until the device or com- ponent has become acclimatised.</li> </ul>
	<ul> <li>Prevent moisture from penetrating into sensitive components.</li> </ul>
	<ul> <li>If moisture forms, switch off the device immediately and allow it to dry properly.</li> </ul>
	Before transporting, fasten the transport lock and disconnect the device from the mains socket.
	The transport temperature must be between -20 °C and +60 °C.
	Humidity must not be condensing. Humidity must be between 10% and 80%.
	Be aware of the weight of the device.
	When transporting using a transport aid (e.g., a pallet truck), the transport aid must be able to carry at least 1.6 times the transport weight of the device.
	Secure the device to prevent it tipping over and falling down during transport.
	Never transport the device sideways or upside down.
Storage conditions	The device must be stored in the original packaging.
	Only store the device in dry rooms.
	The storage temperature must be between -20 °C and +60 °C.
	Humidity must not be condensing. Humidity must be between 10% and 80%.

# 4.2 Transport dimensions and transport weight

Article	Packaging dimensions in mm (L x W x H)	Weight in kg
MIKRO 2.0 / MIKRO 2.0 R	240 x 389 x 255 / 240 x 538 x 250	approx. 20 kg / approx. 33 kg

# 4.3 Fastening the transport lock

## Personnel:

- Trained user
- **1.** Open the lid.

- **2.** Insert the foam cylinder.
- **3.** Close the lid.

## 5 Commissioning

5.1 Unpacking the centrifuge

## CAUTION

Danger of crushing due to parts falling out of the transport packaging.

- Keep the device balanced during the unpacking process.
- Only open the packaging at the points provided for this purpose.

#### CAUTION

Risk of injury from lifting heavy loads.

- Provide an adequate number of helpers.
- Note the weight. See .

# NOTICE

#### Damage to the device due to improper lifting.

- Do not lift the centrifuge by the control panel or the control panel holder.

## 5.2 Removing the transport lock

#### Personnel:

- Trained user
- **1.** Open the lid.
- **2.** Remove the foam cylinder.
- **3.** Store the foam cylinder in a safe place.
- **4.** Close the lid.

## 5.3 Setting up and connecting the centrifuge

Setting up the centrifuge

#### **WARNING**

Operation of the unit in an explosive, corrosive or radioactively contaminated atmosphere is not in accordance with the intended use.

# WARNING

#### **Risk of injury**

Due to insufficient clearance from the centrifuge.

- As per EN / IEC 61010-2-020, no persons, hazardous materials or objects may be present within a safety zone of 300 mm around the centrifuge during a centrifugation run.
- Maintain a distance of 300 mm from the ventilation slots and ventilation openings of the centrifuge.
- The centrifuge's ventilation openings must never be blocked.

## Risk of crushing and damage

Changes in position caused by vibration may cause the device to fall.

- Place the device on a stable and level surface.
- Select an installation area that is capable of bearing the weight of the device.
- Observe national and local regulations on safety and accident prevention.

#### NOTICE

# Damage

Any deviation from the external temperature conditions will result in damage to the samples and device.

- Comply with the maximum and minimum permissible ambient temperatures.
- Avoid placing the device next to heat sources.
- Do not place the device in direct sunlight.
- Protect the device from frost.
- Maintain the required clearance around the device.

#### Personnel:

- Trained user
- **1.** Place the device on a stable and level surface.
- 2. Maintain a distance of 300 mm around the device.
- 3. Comply with the ambient conditions in the technical data ().

#### Connecting the centrifuge

## NOTICE

#### Damage

Property damage caused by unauthorised personnel.

- Do not allow persons without the appropriate authorisation to perform any work on or make changes to devices.
- Only authorised personnel may carry out maintenance and repairs.
- Obtain the manufacturer's approval or guidance before carrying out any work on the device.

## NOTICE

Danger due to condensation in the event of temperature differences

Moisture may damage electrical components.

- Ensure that all surfaces are dry prior to commissioning or maintenance.
- If the temperature changes, wait until the device or component has become acclimatised.
- Prevent moisture from penetrating into sensitive components.
- If moisture forms, switch off the device immediately and allow it to dry properly.

#### Personnel:

- Trained user
- 1. A type B residual current circuit breaker must be used if the device is additionally protected with a residual current circuit breaker in the building installation.

When using a different type, the residual current circuit breaker may either not switch off the unit if there is a fault on the unit, or it may switch off the unit even though there is no fault on the unit.

- 2. Check whether the mains voltage and mains frequency match the specification on the rating plate.
- **3.** Connect the device to a standard mains socket using the mains cable.

## 5.4 Switching on for the first time

#### Personnel:

- Authorisation level 3
- 1. Set the mains switch to [/].
  - ➡ The company logo is briefly displayed.

Animation and introductory text are briefly displayed.

*'Initial start-up'* is displayed.

'Start configuration' is highlighted in blue.

- 2. Press [Rotary, push button].
  - ➡ Language selection is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired language is highlighted in blue.
- 4. Press [Rotary, push button].
  - ➡ Date selection is displayed.
- 5. Rotate [Rotary, push button] until the desired date format setting is selected.
- 6. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.
- 7. Rotate *[Rotary, push button]* until the desired value for the day is selected.
- 8. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.

- 9. Rotate *[Rotary, push button]* until the desired value for the month is selected.
- 10. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.
- **11.** Rotate *[Rotary, push button]* until the desired value for the year is selected.
- **12.** Press [Rotary, push button].
  - Date selection is displayed.
- **13.** Rotate *[Rotary, push button]* until the desired time format setting is selected.
- 14. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.
- **15.** Rotate *[Rotary, push button]* until the desired value for the hours is selected.
- **16.** Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.
- 17. Rotate [*Rotary, push button*] until the desired value for the minutes is selected.
- 18. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.
- **19.** Rotate *[Rotary, push button]* until the desired value for the seconds is selected.
- 20. Press [Rotary, push button].
  - ➡ PIN entry is displayed.
- **21.** Rotate *[Rotary, push button]* and confirm by pressing to select the desired number. Repeat the process until the desired 8 numbers have been selected.
  - ➡ ✓ symbol is highlighted in blue.
- 22. Press [Rotary, push button].
  - ➡ The settings are stored.

'Initial start-up successful' is displayed.

'Retart centrifuge' is highlighted in blue.

- **23.** Press [Rotary, push button].
  - Centrifuge is restarted.

## 5.5 Switching the centrifuge on and off.

Switching the centrifuge on Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

22 / 77

- ▶ Set the mains switch to [/].
  - ➡ The following displays appear one after the other:
    - Progress bar
    - Company logo
  - ➡ Quick Adjust menu is selected.

	$\bigcirc$
۵	
0	

In the Settings menu, the user can choose whether to always start in the Program menu or the Quick Adjust menu. You can also choose whether to start with the last menu that was executed. See .

To simplify and clarify presentation, in this document it is always assumed that the Quick Adjust menu is started.

#### Switching off the centrifuge

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The rotor is stationary.

- \_\_\_ Set the mains switch to [0].
  - Device shuts down.

Blue background with symbol is displayed.

## 6 Operation

#### WARNING

Flammable gas (R290) - possible leakage from the centrifuge

Failure to observe this can lead to a flash fire, which can result in minor to moderate injury and damage.

- No ignition sources or naked flames in the room.

## 6.1 Opening and closing the lid

Opening the lid

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The centrifuge is switched on.

The rotor is stationary.

- Press the *[STOP, QUICK STOP, OPEN LID]* button.
  - ➡ The lid unlocks by means of a motor.

Closing the lid

# AUTION

#### Risk of crushing

Danger of crushing fingers when the closing motor presses the lid against the seal.

- Keep body parts clear of the lid danger zone when closing the lid.
- Press the lid down from above to close it.

## NOTICE

## Damage

Slamming the lid damages the appliance.

- Close the lid slowly.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
  - Close the lid and press the front edge of the lid down gently.
    - ➡ The lid locks using a motor.

## 6.2 Removing and inserting the rotor

#### Rotor removal

#### Personnel:

- Trained user
- 1. Den the centrifuge lid.
- **2.** Grip beneath the removal aid (1) with your index finger and middle finger.
- **3.** Press the release pin (2) down with your thumb.
- **4.** Pull the removal aid (1) upwards.
- 5. Lift the rotor (5) with the rotor lid (4) upwards from the rotor mount (6) and remove.

1	Removal aid	4	Rotor lid
2	Release pin	5	Rotor
3	Lid release	6	Rotor mount

Inserting the rotor

#### Personnel:

Trained user

The centrifuge lid is open.

- 1. Clean the rotor mount (6) and the inner surface of the rotor adapter (XXX) for the rotor.
- **2.** Grip beneath the removal aid (1) with your index finger and middle finger.
  - ➡ Removal aid (xxxx) moves upwards.
- **3.** Place the rotor (5) vertically on the rotor mount (6).
- **4.** Check whether the rotor (5) is correctly inserted.

## 6.3 Fitting and removing the rotor lid

Removing the rotor lid

## Personnel:

Trained user

The lid is open.

**1.** Press the lid release (3) together with your index finger and thumb.

2. Lift the rotor lid (4) upwards from the rotor (5) and remove.

Fig. 5: Fitting and removing the rotor lid

1	Removal aid	4	Rotor lid
2	Release pin	5	Rotor
3	Lid release	6	Rotor mount

#### Fitting the rotor lid

#### Personnel:

Trained user

The lid is open.

- **1.** Press the lid release (3) together with your index finger and thumb.
- 2. Place the rotor lid (4) vertically on the rotor (5).
- **3.** Check whether the rotor lid (4) is correctly engaged.

Check whether the lid release (3) is in the correct position.

## 6.4 Loading

Filling centrifuge tubes

#### WARNING

#### Risk of injury from contaminated sample material

Leakage of contaminated sample material from the sample tube during centrifugation.

- When working with hazardous substances, only use centrifuge tubes with special screw caps.
- For materials in risk groups 3 and 4, additionally use a biosafety system in accordance with the WHO 'Laboratory Biosafety Manual' alongside sealable centrifuge tubes.

## NOTICE

#### Damage

Device damage due to corrosive substances.

- Do not centrifuge highly corrosive substances.
- Do not allow corrosive substances to enter the device.
- Check the device regularly for corrosion.

Standard glass centrifuge tubes can be loaded up to RCF 4000 (DIN 58970 part 2).

#### Personnel:

- Trained user
- Fill centrifuge tubes outside the centrifuge.

The maximum capacity of the centrifuge tubes specified by the manufacturer must not be exceeded.

With angle rotors, the centrifuge tubes must only be filled to the extent that no liquid can be ejected from the tubes during the centrifugation run.

It must be ensured that there is a uniform fill level in the tubes in order to keep the weight differences in the centrifuge tubes as low as possible.

**Symbols on the angle rotor** The following symbols appear on the rotor to help with even rotor loading:

Fig. 6: Rotor with symbols

- 1 Division of the rotor into four segments
- 2 Division of the rotor into three segments

# Loading the angle rotors

- Division of the rotor into four segments. Two thick arcs are used for marking. Loading aid for even distribution of the centrifugation tubes over four rotor segments.
- Division of the rotor into three segments.
   Location numbers are used for marking, and these appear in a larger font.

Loading aid for even distribution of the centrifugation tubes over three rotor segments.

# Personnel:

- Trained user
- 1. Check that the rotor is firmly seated.
- **2.** The centrifuge tubes must be distributed evenly over all locations on the rotor.

No liquid must be allowed to enter the rotor and the centrifuging chamber when loading the rotor.

With rotors, the centrifuge tubes must only be filled to the extent that no liquid can be ejected from the tubes during the centrifugation run.

The weight of the permissible filling capacity is indicated on each rotor. The weight must not be exceeded.

# 6.5 Centrifugation parameters

# 6.5.1 Relative centrifugal force, RCF

The relative centrifugal force (RCF) is dependent on the speed and the centrifuging radius.

RCF is specified as a multiple of the acceleration due to gravity (g).

The RCF is a dimensionless numerical value and is used to compare the separation and sedimentation performance.

$$\mathrm{RCF} = \left(\frac{\mathrm{RPM}}{1000}\right)^2 * \mathrm{r} * 1,118$$

$$RPM = \sqrt{\frac{RCF}{r * 1,118}} * 1000$$

RCF = Relative Centrifugal Force

RPM = revolutions per minute,  $[RPM] = min^{-1}$ .

r = distance from the axis of rotation to the bottom of the centrifuge tube, [r] = mm.

# 6.5.2 Centrifugation of substances or mixtures of substances with a density higher than 1.2 kg/dm<sup>3</sup>

The density of the substances or mixtures of substances must not exceed 1.2 kg/dm<sup>3</sup> during centrifugation at maximum speed. The speed must be reduced for substances or substance mixtures with a higher density. The permissible speed can be calculated using the following formula:

Reduced speed  $(n_{red}) = \sqrt{\frac{1,2}{\text{Greater density [kg/dm<sup>3</sup>]}}} * \text{maximum speed [RPM]}$ 

For example: Maximum speed 4000 RPM, density 1.6 kg/dm<sup>3</sup>

$$n_{red} = \sqrt{\frac{1,2(\text{kg/dm}^3)}{1,6(\text{kg/dm}^3)}} * 4000 \text{ RPM} = 3464 \text{ RPM}$$

If, in exceptional cases, the maximum load indicated on the bucket is exceeded, the speed must also be reduced. The permissible speed can be calculated using the following formula:

Reduced speed  $(n_{red}) = \sqrt{\frac{\text{maximum load [g]}}{\text{actual load [g]}} * \text{maximum speed [RPM]}}$ 

For example: Maximum speed 4000 RPM, maximum load 300 g, actual load 350 g

$$n_{red} = \sqrt{\frac{300 \text{ g}}{350 \text{ g}}} * 4000 \text{ RPM} = 3703 \text{ RPM}$$

Please contact the manufacturer if you are not sure.

3.5.3 Ramp-up and brake levels				
		1		
		2 3 4		
		1 2 3 4 5 6 7 8 9 10 10 9 8 7 6 5 4 3 2 1 0 9 8 7 6 5 4 3 2 1 0 9 8 7 6 5 4 3 2 1		
		<ul> <li>Fig. 7: Ramp-up and brake levels</li> <li>Runtime (runtime measurement after the start of the centrifugation run)</li> <li>Ramp-up levels</li> <li>Runtime (runtime measurement after reaching the speed)</li> <li>Brake levels</li> </ul>		
Runtime		You can set how the runtime is measured:		
		Runtime measurement starts after the start of the centrifugation run (1).		
		Runtime measurement starts after reaching centrifugation speed $(\mathcal{3})$ .		
		To set the runtime measurement, see or .		
Ramp-up	evels	10 ramp-up levels can be set (2).		
		Ramp-up level 1: longest time until the speed is reached.		
		Ramp-up level 10: shortest time until the speed is reached.		
		To set the ramp-up levels, see .		
Brake levels		11 brake levels can be set $(4)$ .		
		Brake level 10: shortest time until the rotor comes to a standstill.		
		Brake level 1: second longest time until the rotor comes to a standstill.		
		Brake level 0: unbraked ramp-down of the rotor. Longest time until the rotor comes to a standstill.		
		To set the brake levels, see .		
0.0				
6.6 U	iser interface			
6.6.1	Dealing with message	jes		
		An information, warning or error message is displayed in the information line of the user interface if the centrifuge is in a state that the user needs to be informed about.		

For certain messages, both an audible signal (see ) and a visual signal (see ) can also be set.

Warning messages cannot be acknowledged by level 1 users.

### Information about the centrifugation run

12345123451234512345

- Show details

	1. Press the [INFO] button.
	Details of the information about the centrifugation run are displayed on the user interface.
	2. Press the [BACK] button.
	Details of the information about the centrifugation run are hidden.
	<b>Option</b> Remove message and information in the detailed view by pressing <i>[Rotary, push button]</i> or opening the lid.
Warning messages	12345123451234512345 123451234512345
- Show details	Warning message appears in the information line.
	<ul> <li>Details of the warning message are displayed on the user interface</li> </ul>
	2. ► Press the <i>/BACK</i> / button.
	<ul> <li>Details of the warning message are hidden.</li> </ul>
- Acknowledge message	Warning message appears in the information line.
	1. [Rotary, push button] or press the [BACK] button.
	Query is displayed.
	2. Press [Rotary, push button].
	Message is acknowledged.
	A PIN must be entered for authorisation level 1.
	It may be necessary to enter a PIN for authorisation level 2.
Error messages	A 123451234512345123451
	- 2345123451234512345
- Show details	Error message appears in the information line.
- Show details	Error message appears in the information line.  1. Press the <i>[INFO]</i> button.
- Show details	2345123451234512345         Error message appears in the information line.         1. Press the [INFO] button.            → Details of the error message are displayed on the user interface.
- Show details	<ul> <li>Error message appears in the information line.</li> <li>1. Press the <i>[INFO]</i> button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>2. Press the <i>[BACK]</i> button.</li> </ul>
- Show details	<ul> <li>Error message appears in the information line.</li> <li>1. Press the [/NFO] button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>2. Press the [BACK] button.</li> <li>Details of the error message are hidden.</li> </ul>
- Show details - Acknowledge message	<ul> <li>Error message appears in the information line.</li> <li>1. Press the <i>[INFO]</i> button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>2. Press the <i>[BACK]</i> button.</li> <li>Details of the error message are hidden.</li> <li>Error message appears in the information line.</li> </ul>
- Show details - Acknowledge message	<ul> <li>Error message appears in the information line.</li> <li>1. Press the [/NFO] button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>2. Press the [BACK] button.</li> <li>Details of the error message are hidden.</li> <li>Error message appears in the information line.</li> <li>Switching the centrifuge off and on again, see - Chapter 5.5</li> </ul>
- Show details - Acknowledge message	<ul> <li>Error message appears in the information line.</li> <li>1. Press the [/NFO] button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>2. Press the [BACK] button.</li> <li>Details of the error message are hidden.</li> <li>Error message appears in the information line.</li> <li>Switching the centrifuge off and on again, see - Chapter 5.5 'Switching the centrifuge on and off.' on page 22.</li> <li>If the error message persists, contact the manufacturer's service.</li> </ul>
- Show details - Acknowledge message	<ul> <li>Error message appears in the information line.</li> <li>1. Press the <i>[INFO]</i> button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>2. Press the <i>[BACK]</i> button.</li> <li>Details of the error message are hidden.</li> <li>Error message appears in the information line.</li> <li>Switching the centrifuge off and on again, see - <i>Chapter 5.5 'Switching the centrifuge on and off.' on page 22.</i></li> <li>If the error message persists, contact the manufacturer's service department.</li> </ul>
- Show details - Acknowledge message	<ul> <li>Error message appears in the information line.</li> <li>1. Press the <i>[INFO]</i> button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>2. Press the <i>[BACK]</i> button.</li> <li>Details of the error message are hidden.</li> <li>Error message appears in the information line.</li> <li>Switching the centrifuge off and on again, see - <i>Chapter 5.5 'Switching the centrifuge on and off.' on page 22.</i></li> <li>If the error message persists, contact the manufacturer's service department.</li> </ul>
<ul> <li>Show details</li> <li>Acknowledge message</li> <li>6.6.2 Basic displays</li> </ul>	<ul> <li>Error message appears in the information line.</li> <li>1. Press the [INFO] button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>2. Press the [BACK] button.</li> <li>Details of the error message are hidden.</li> <li>Error message appears in the information line.</li> <li>Switching the centrifuge off and on again, see - Chapter 5.5 'Switching the centrifuge on and off.' on page 22.</li> <li>If the error message persists, contact the manufacturer's service department.</li> </ul>
<ul> <li>Show details</li> <li>Acknowledge message</li> <li>6.6.2 Basic displays</li> </ul>	<ul> <li>Error message appears in the information line.</li> <li>1. Press the <i>[INFO]</i> button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>2. Press the <i>[BACK]</i> button.</li> <li>Details of the error message are hidden.</li> <li>Error message appears in the information line.</li> <li>Switching the centrifuge off and on again, see - Chapter 5.5 <i>'Switching the centrifuge on and off.' on page 22.</i></li> <li>If the error message persists, contact the manufacturer's service department.</li> </ul>
<ul> <li>Show details</li> <li>Acknowledge message</li> <li>6.6.2 Basic displays</li> </ul>	<ul> <li>Error message appears in the information line.</li> <li>Press the <i>[INFO]</i> button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>Press the <i>[BACK]</i> button.</li> <li>Details of the error message are hidden.</li> <li>Error message appears in the information line.</li> <li>Switching the centrifuge off and on again, see <i>Chapter 5.5 'Switching the centrifuge on and off.' on page 22.</i></li> <li>If the error message persists, contact the manufacturer's service department.</li> </ul>
<ul> <li>Show details</li> <li>Acknowledge message</li> <li>6.6.2 Basic displays</li> </ul>	<ul> <li>Error message appears in the information line.</li> <li>Press the <i>[INFO]</i> button.</li> <li>Details of the error message are displayed on the user interface.</li> <li>Press the <i>[BACK]</i> button.</li> <li>Details of the error message are hidden.</li> <li>Error message appears in the information line.</li> <li>Switching the centrifuge off and on again, see - Chapter 5.5 <i>Switching the centrifuge on and off.' on page 22.</i></li> <li>If the error message persists, contact the manufacturer's service department.</li> </ul>

Scroll bar

position of the selected entry.

In the various menus, the scroll bars indicate the length of the list and the

# Status bar

The following symbols can be displayed in the status bar:

Symbol	Meaning
•	Indicates the progress of the centrifugation run.
	Start of the centrifugation run as soon as the lid is closed.
	For uncooled centrifuges only: Lid opens automatically after a successful centrifuga- tion run
\$3	Authorisation level Level 1: Restricted rights Level 2: Standard rights Level 3: Administrator
!	The set parameters do not match the rotor used.

The following symbols can be displayed in the information lines:

Symbol	Meaning
$\overline{\Box}$	The lid is closed.
$\Box$	The lid is open.
í	More information can be displayed by pressing the <i>[INFO]</i> button.
$\triangle$	error message
X	Wait
	Short-term centrifugation
	Сору
ক্ট্য	Settings

Symbol	Meaning
~	Action performed successfully.
×	Action cancelled.

#### 6.6.3 Keypad

	$ \begin{array}{cccccccccccccccccccccccccccccccccccc$
	<ul> <li>Fig. 8: Keypad</li> <li>1 Country-specific keypad</li> <li>2 Delete input</li> <li>3 Space bar</li> <li>4 Toggle between upper and lower case letters</li> <li>5 Special characters</li> <li>6 Letters and numbers</li> </ul>
Letters and numbers	<ol> <li>Rotate [Rotary, push button] until the desired letter or number is high-lighted in blue.</li> <li>Press [Rotary, push button].</li> <li>The letter or number is transferred to the input field.</li> </ol>
Space bar	<ol> <li>Rotate [Rotary, push button] until the space bar (3) is highlighted in blue.</li> <li>Press [Rotary, push button].</li> </ol>

- ➡ A space is transferred to the input field.
- **1.** Rotate *[Rotary, push button]* until the symbol () is highlighted in blue.
- 2. Press [Rotary, push button].
  - ➡ Upper or lower case letters are displayed.
- **1.** Rotate *[Rotary, push button]* until the symbol () is highlighted in blue.
- 2. Press [Rotary, push button].
  - ➡ The special characters are displayed.
- 3. Rotate [Rotary, push button] until the desired special character is highlighted in blue.

Toggle between upper and

lower case letters

**Special characters** 

	<ul> <li>4. Press [Rotary, push button].</li> <li>→ The special character is transferred to the input field.</li> </ul>
Delete input	<ol> <li>Rotate [Rotary, push button] until the symbol () is highlighted in blue.</li> <li>Press [Rotary, push button].</li> <li>The last entry is deleted from the input field.</li> </ol>
Country-specific keypad	<b>1.</b> Rotate <i>[Rotary, push button]</i> until the language code of the country () is highlighted in blue.
	<ul> <li>Press [Rotary, push button].</li> <li>♦ The country-specific keypad for the displayed country is displayed.</li> </ul>

 The next possible language code () is displayed in alphabetical order.

# 6.7 Menus

# 6.7.1 Main menu

# Fig. 9: Main menu

- 1 Quick Adjust menu
- 2 Program menu
- 3 Settings menu
- 4 Status bar

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- Press [Rotary, push button].
  - ➡ The background of the selected menu symbol turns white.
    - Menu contents are displayed.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

A menu is selected.

- **1.** Press the *[BACK]* button.
  - The background of the menu symbol turns blue.

Menu contents are hidden.

Selecting a menu

Changing the menu

- 2. Rotate [Rotary, push button].
  - ➡ The background of the next menu symbol turns blue.
- 3. Press [Rotary, push button].
  - The background of the next menu symbol turns white.
     Menu contents are displayed.

# 6.7.2 Quick Adjust menu

Fig. 10: Quick Adjust menu

- 1 Centrifuging radius, speed (RPM) or relative centrifugal force (RCF)
- 2 Runtime
- 3 Temperature (only for devices with cooling)
- 4 Ramp-up and brake level

#### Selection of parameters

- Press [Rotary, push button].
  - ➡ The background of the next menu symbol turns blue.

# 6.7.3 Program menu

#### Fig. 11: Program menu

- 1 Number of programs created / maximum number of possible programmes
- 2 Create new program

1. Rotate [Rotary, push button].

### Selection of programs

➡ Navigate to the desired program.

- **2.** Press [Rotary, push button].
  - ➡ Program data is displayed.

6.7.4 Settings menu

Selection of submenus

- Fig. 12: Settings menu
- 1. Rotate [Rotary, push button].
   Navigate to the desired menu item.
- 2. Press [Rotary, push button].
  - ➡ Menu is selected.
- 6.8 Centrifugation
- 6.8.1 Continuous operation

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The rotor is installed and the centrifuge lid is closed.

In the Quick Adjust menu or in the program menu, the symbol  $\ensuremath{\bigcirc}\xspace$  is set during runtime.

- 1. Press the *[START]* button.
  - ➡ The centrifugation run is started.

The following summary is displayed during the centrifugation run:

(1)		
2	- (?) mm RCF 999 99 999	mm RCF 999   99 999
3—	- Ō 99:59:59	∕-   <sup>h:m:s</sup> 99:59:59
4	€ °C 999	°⊂ 999
5	$\neg$	99 99
(6)		【靣 台 \$1

- 1 Selected program
- 2 Current centrifuging radius, speed (RPM) or relative centrifugal force (RCF)
- 3 Current runtime
- 4 Current temperature (for devices with cooling)
- 5 Pre-set ramp-up level and brake level
- 6 Status bar
- 2. Press the [STOP, QUICK STOP, OPEN LID] button to end the centrifugation run.
  - ➡ Ramp-down takes place with the set brake level.

The lid can be opened shortly after the rotor has come to a standstill.

Information line with information about the centrifugation run is displayed.

- 3. If required: Press the *[INFO]* button after the centrifugation run.
  - Centrifugation run parameters are displayed.
- **4.** If required: Press the *[BACK]* button.
  - Centrifugation run parameters are hidden.
- 5. Opening the lid
  - or

Press the [BACK] button.

➡ The original screen is displayed.

# 6.8.2 Centrifugation with time preselection

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The rotor is installed and the centrifuge lid is closed.

The desired parameters are set in the Quick Adjust menu or the program menu.

- 1. Press the *[START]* button.
  - ➡ The centrifugation run is started.

The following summary is displayed during the centrifugation run:

(1)		
ğ—		mm RCF 999   99 999
3—	Ō h:m:s 99:59:59	✓   <sup>h:m:s</sup> 99:59:59
4	°C 999	° <b>c</b> 999
5		99 99
6		靣 ՝ े \$1

- 1 Selected program
- 2 Current centrifuging radius, speed (RPM) or relative centrifugal force (RCF)
- 3 Current runtime
- 4 Current temperature (for devices with cooling)
- 5 Pre-set ramp-up level and brake level
- 6 Status bar
- 2. Press the *[STOP, QUICK STOP, OPEN LID]* button to cancel the centrifugation run.
  - ➡ Ramp-down takes place with the set brake level.

The lid can be opened shortly after the rotor has come to a standstill.

Information line with information about the centrifugation run is displayed.

- **3.** If required: Press the *[INFO]* button after the centrifugation run.
  - Centrifugation run parameters are displayed.
- **4.** If required: Press the *[BACK]* button.
  - ➡ Centrifugation run parameters are hidden.
- 5. Opening the lid

or

Press the *BACK* button.

➡ The original screen is displayed.

# 6.8.3 Short-term centrifugation

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The rotor is installed and the centrifuge lid is closed.

The desired parameters are set in the Quick Adjust menu.

- 1. Press and hold the [SHORT-TERM CENTRIFUGATION] button.
  - ➡ The centrifugation run is started.

Short-term centrifugation can be started when the Quick Adjust menu, the main menu or the program list is displayed.

The following summary is displayed during the centrifugation run:

- 1 Current temperature (for devices with cooling)
- 2 Current runtime
- 3 Current centrifuging radius, speed (RPM) or relative centrifugal force (RCF)
- 4 Short-term centrifugation indicator
- 5 Pre-set ramp-up level
- 6 Pre-set speed (RPM) or relative centrifugal force (RCF)
- 7 Pre-set brake level
- 2. Release the *[SHORT-TERM CENTRIFUGATION]* button to end the centrifugation run.
  - Ramp-down takes place with the set brake level.

The lid can be opened shortly after the rotor has come to a standstill.

Information line with information about the centrifugation run is displayed.

- 3. If required: Press the [INFO] button.
  - Centrifugation run parameters are displayed.
- 4. If required: Press the [BACK] button.
  - ➡ Centrifugation run parameters are hidden.
- 5. Opening the lid
  - or
  - Press the [BACK] button.
  - ➡ The original screen is displayed.

# 6.9 Quick stop function

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

▶ Press the *[STOP, QUICK STOP, OPEN LID]* button twice

Press and hold the [STOP, QUICK STOP, OPEN LID] button.

Ramp-down takes place at brake level "10" (shortest ramp-down time).

# 6.10 Software operation

6.10.1 Set Quick Adjust

Quick Adjust menu is selected.

### Personnel:

or

- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* to change the value.
- 2. Press [Rotary, push button].
  - ➡ Value is saved.

The background of the next menu symbol turns blue.

The runtime is displayed in bold and can be changed.

3. Rotate [Rotary, push button] to change the value.

For centrifuging in continuous operation, set the  $\bigcirc$  symbol. The symbol is located between *'99:59:59 h:m:s'* and *'00:00:01 h:m:s'*.

- 4. Press [Rotary, push button].
  - ➡ Value is saved.

The background of the next menu symbol turns blue.

For centrifuges with cooling: The temperature is displayed in bold and can be changed.

For centrifuges without cooling: The ramp-up level is displayed in bold and can be changed. The next two steps are skipped.

- 5. Rotate [Rotary, push button] to change the value.
- 6. Press [Rotary, push button].
  - ➡ Value is saved.

The background of the symbol for the ramp-up and brake level turns blue.

The ramp-up level is displayed in bold and can be changed.

- 7. Rotate [Rotary, push button] to change the value.
- 8. Press [Rotary, push button].
  - ➡ The brake level is displayed in bold and can be changed.
- 9. Rotate [Rotary, push button] to change the value.
- **10.** Press the *[BACK]* button.

➡ The changes are stored.

#### Changing the relative centrifugal force (RCF), centrifuging

- Authorisation level 2
- Authorisation level 3

Speed, relative centrifugal force (RCF), centrifuging radius are selected.

radius

- 1. Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
- 2. Rotate [Rotary, push button] to change the value.
  - ➡ Selected value is displayed in bold.
- 3. Press [Rotary, push button].
  - ➡ Radius is selected and displayed in bold.
- **4.** Rotate [*Rotary, push button*] to change the value.
- 5. Press the [BACK] button.
  - ➡ The changes are stored.

# Personnel:

- Authorisation level 2
- Authorisation level 3
- Runtime is selected.
- 1. [CONTEXT BUTTON, TOGGLE BUTTON] button
- 2. Rotate [Rotary, push button] to change the value.
  - ➡ Selected value is displayed in bold.
- **3.** Press the *[BACK]* button.
  - ➡ The changes are stored.

Changing parameters during the centrifugation run

#### Personnel:

- Authorisation level 2
- Authorisation level 3

Centrifugation run in the Quick Adjust menu is active.

The runtime display is set to 'classic'.

- 1. Press and hold the *[Rotary, push button]* button for 5 seconds.
  - The speed (RPM) or relative centrifugal force (RCF) parameters of the current centrifugation run are shown in bold in the Quick Adjust menu.
- **2.** Rotate [*Rotary, push button*] to change the value.
- **3.** Press [Rotary, push button].
  - The runtime parameters of the current centrifugation run are shown in bold in the Quick Adjust menu.
- **4.** Rotate *[Rotary, push button]* to change the value.
- 5. Press [START] to save the changes to the parameters in Quick Adjust.

Press [BACK] to discard the changes to the parameters in Quick Adjust.

Displaying parameters of the last centrifugation run

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- Press the *[INFO]* button.

Rev. -- / 05.2025

The parameters of the last centrifugation run are displayed in the Quick Adjust menu, the program list (only when the program names are displayed) and the main menu.

# Changing the runtime

# 6.10.2 Program menu for authorisation levels

Creating a program

#### Personnel:

Authorisation level 3

Program menu is selected.

- **1.** Rotate *[Rotary, push button]* to the right until the create new program symbol is highlighted in blue.
- **2.** Press [Rotary, push button].
  - ➡ Keyboard opens.
- **3.** Rotate and press *[Rotary, push button]* to enter the program name.
- 4. Press the [BACK] button.
  - Program is saved with the parameters from the Quick Adjust menu.

### Retrieving a program

#### Personnel:

- Authorisation level 2
- Authorisation level 3

Program menu is selected.

- **1.** Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
- 2. Press [Rotary, push button].
  - ➡ Program parameters are displayed.
- 3. Press [Rotary, push button].
  - ➡ Program parameters are opened.

#### Changing a program

Program menu is selected.

Program is called.

Fig. 13: Program mode menu - Change program

1 Centrifuging radius, speed (RPM) or relative centrifugal force (RCF)

- 2 Runtime
- 3 Temperature (only for devices with cooling)
- 4 Ramp-up and brake level

#### Personnel:

- Authorisation level 3
- 1. Rotate [Rotary, push button] to change the value.
- 2. Press [Rotary, push button].
  - Value is saved.

The background of the next menu symbol turns blue. The runtime is displayed in bold and can be changed.

3. Rotate [Rotary, push button] to change the value.

For centrifuging in continuous operation, set the  $\bigcirc$  symbol. The symbol is located between *'99:59:59 h:m:s'* and *'00:00:01 h:m:s'*.

- 4. Press [Rotary, push button].
  - Value is saved.

The background of the next menu symbol turns blue.

For centrifuges with cooling: The temperature is displayed in bold and can be changed.

For centrifuges without cooling: The ramp-up level is displayed in bold and can be changed. The next two steps are skipped.

- 5. Rotate [Rotary, push button] to change the value.
- 6. Press [Rotary, push button].
  - ➡ Value is saved.

The background of the symbol for the ramp-up and brake level turns blue.

The ramp-up level is displayed in bold and can be changed.

- 7. Rotate [Rotary, push button] to change the value.
- 8. Press [Rotary, push button].
  - ➡ The brake level is displayed in bold and can be changed.
- 9. Rotate [Rotary, push button] to change the value.
- **10.** Press the *[BACK]* button.
  - ➡ The changes are stored.

#### Personnel:

Authorisation level 3

Speed, relative centrifugal force (RCF), centrifuging radius are selected.

- 1. Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
- 2. Rotate [Rotary, push button] to change the value.
  - Selected value is displayed in bold.
- **3.** Press [*Rotary, push button*].
  - ➡ Radius is selected and displayed in bold.
- **4.** Rotate *[Rotary, push button]* to change the value.
- 5. Press the [BACK] button.
  - ➡ The changes are stored.

#### Personnel:

Authorisation level 3

Runtime is selected.

1. [CONTEXT BUTTON, TOGGLE BUTTON] button

Changing speed, relative centrifugal force (RCF), centrifuging radius

Changing the runtime

- 2. Rotate [Rotary, push button] to change the value.
  - Selected value is displayed in bold.
- **3.** Press the *[BACK]* button.
  - ➡ The changes are stored.

#### Creating program favourites

#### Personnel:

Authorisation level 3

Program menu is selected.

- **1.** Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
- 2. Press [Rotary, push button].
  - ➡ Program parameters are displayed.
- 3. Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
  - ➡ Submenu appears.
- **4.** Rotate *[Rotary, push button]* until the Favourites submenu symbol is highlighted in blue.
- 5. Press [Rotary, push button].
  - ➡ The Favourites submenu is displayed.

Favourite places that are already assigned are marked with a solid star.

- 6. Rotate [Rotary, push button] until the desired free favourite place is highlighted in blue.
- 7. Press [Rotary, push button].
  - ➡ Program is assigned a favourite place.

#### Deleting program favourites

#### Personnel:

Authorisation level 3

Program menu is selected.

- **1.** Rotate *[Rotary, push button]* until the desired favourite program is highlighted in blue.
- 2. Press [Rotary, push button].
  - Program parameters are displayed.
- 3. Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
  - ➡ Submenu appears.
- **4.** Rotate *[Rotary, push button]* until the Favourites submenu symbol is highlighted in blue.
- 5. Press [Rotary, push button].
  - ➡ The Favourites submenu is displayed.

The assigned favourite place is highlighted in blue.

- 6. Press [Rotary, push button].
  - Program is removed from favourite place.

Copying a program

#### Personnel:

Authorisation level 3

Program menu is selected.

- **1.** Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
- 2. Press [Rotary, push button].
  - ➡ Program parameters are displayed.
- 3. Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
  - Submenu appears.
- **4.** Rotate *[Rotary, push button]* until the Copy submenu symbol is highlighted in blue.
- 5. Press [Rotary, push button].
  - Keyboard opens.
- 6. Rotate and press *[Rotary, push button]* to enter the program name.
- 7. Press the [BACK] button.
  - Program is saved with the copied parameters.
     Information line indicates successful execution of the action.

Renaming a program

#### Personnel:

Authorisation level 3

Program menu is selected.

- **1.** Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
- **2.** Press [Rotary, push button].
  - Program parameters are displayed.
- 3. Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
  - Submenu appears.
- **4.** Rotate *[Rotary, push button]* until the Rename submenu symbol is highlighted in blue.
- 5. Press [Rotary, push button].
  - Keyboard opens.
- 6. Rotate and press [Rotary, push button] to enter the program name.
- 7. Press the [BACK] button.
  - ➡ Program is saved with the new name.

#### Deleting a program

#### Personnel:

Authorisation level 3

Program menu is selected.

- **1.** Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
- **2.** Press [*Rotary, push button*].
  - Program parameters are displayed.
- 3. Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
  - Submenu appears.
- **4.** Rotate *[Rotary, push button]* until the Delete submenu symbol is highlighted in blue.
- 5. Press [Rotary, push button].
  - Confirmation query appears
- 6. Rotate and press [Rotary, push button] to confirm the query.

7. Press [Rotary, push button].

Program is deleted.

# 6.10.3 Device information

System information

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'Device information' submenu is selected in the settings menu.

- **1.** Rotate *[Rotary, push button]* until the *'System information'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.

The following information can be found in the menu:

- Machine name
- Sales number
- Equipment number
- Serial number
- HBI software version
- CPU serial number
- ZS software version
- Manufacturer URL
- Next service date (date/time (UTC))

#### Rotor info

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'Device information' submenu is selected in the settings menu.

- **1.** Rotate *[Rotary, push button]* until the *'Rotor info'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired rotor is highlighted in blue.

- 4. Press [Rotary, push button].
  - Menu is displayed.

The following information can be found in the menu:

- Sales number
- Serial number
- Expiry date
- Machine-dependent, maximum speed
- Maximum rotor cycles, maximum number of rotor cycles
- Current rotor cycles
- Remaining rotor cycles
- Lowest achievable temperature at maximum speed (only for centrifuges with cooling).

Error history

#### Personnel:

- Authorisation level 2
- Authorisation level 3

The 'Device information' submenu is selected in the settings menu.

- **1.** Rotate *[Rotary, push button]* until the *'Error history'* menu is highlighted in blue.
- **2.** Press [Rotary, push button].
  - ➡ Menu is displayed.

Warnings and errors that have occurred are displayed:

- Date of the warning or error
- Time of the warning or error
- Number of the warning or error
- **3.** Rotate *[Rotary, push button]* until the desired error is highlighted in blue.
- **4.** Press [Rotary, push button].
  - Error short text, error number and complete error text are displayed.

#### **Operating hours**

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'Device information' submenu is selected in the settings menu.

- **1.** Rotate *[Rotary, push button]* until the *[Querying]* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.

The following information can be found in the menu:

- Current runtime
- Operating hours connected to the grid
- Total runtime

#### Data export to USB flash drive

In the menu, system log files and audit log files can be exported to a USB flash drive.

The following file systems can be used for USB flash drives:

■ FAT, FAT32, ext2, ext3, ext4

#### Personnel:

- Authorisation level 2
- Authorisation level 3

USB flash drive is inserted.

The 'Device information' submenu is selected in the settings menu.

- 1. Rotate [Rotary, push button] until the 'Data export to USB flash drive' menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired data for export is highlighted in blue.
- 4. Press [Rotary, push button].
  - Menu is displayed.
- 5. Rotate [*Rotary, push button*] until the desired time period for the export is highlighted in blue.
- 6. Press [Rotary, push button].
  - Menu with export requirements is displayed.
- 7. Press [Rotary, push button].
  - ➡ Export is started.

#### Service contact details

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'Device information' submenu is selected in the settings menu.

- **1.** Rotate *[Rotary, push button]* until the *'Service contact details'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - ➡ Menu is displayed.

The following information can be found in the menu: Service contact details:

- Phone:
- E-mail
- Note (information entered by service)

# 6.11 User settings

Starting point in Quick Adjust

The menu option is selected so that the starting point in Quick Adjust is either the RPM/RCF or the time.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- 1. Rotate [Rotary, push button] until the 'Starting point in Quick Adjust' menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired setting is selected.
- 4. Press the [BACK] button.
  - ➡ The settings are stored.

#### **Runtime display**

The menu settings allow you to choose between a *'enlarged'* or *'classic'* display during centrifugation.

#### 'enlarged':

The enlarged runtime display is shown during a centrifugation run or continuous centrifugation. In enlarged view, no changes can be made to the parameters during a Quick Adjust centrifugation run.

'classic':

The classic runtime display is shown during a centrifugation run or during continuous centrifugation.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- 1. Rotate [Rotary, push button] until the 'Runtime display' menu is highlighted in blue.
- **2.** Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired setting is selected.
- **4.** Press the *[BACK]* button.
  - ➡ The settings are stored.

#### Step size for runtime settings

In the menu, the steps are set to 1 second or 15 seconds when setting the runtime.

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- 1. Rotate [*Rotary, push button*] until the [*Step size for runtime settings*] menu is highlighted in blue.

- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired setting is selected.
- 4. Press the [BACK] button.
  - ➡ The settings are stored.

#### Lid opens after centrifugation

The setting "Lid opens after centrifugation" is only available for Mikro 2.0. The menu is used to set whether the centrifuge lid opens automatically after centrifugation is complete.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Lid opens after centrifugation'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired setting is selected.
- 4. Press the [BACK] button.
  - ➡ The settings are stored.

Lid opens when switched on The menu is used to set whether the centrifuge lid opens automatically when the centrifuge is switched on.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *[Lid opens when switched on]* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- 3. Rotate [Rotary, push button] until the desired setting is selected.
- 4. Press the [BACK] button.
  - ➡ The settings are stored.

#### Automatic start after lid closure

In the menu, centrifugation is set to start automatically once the centrifuge lid has been closed.

Automatic start :

Centrifugation only starts when Quick Adjust or the program list is displayed.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Automatic start after lid closure'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- 3. Rotate [Rotary, push button] until the desired setting is selected.
- 4. Press the [BACK] button.
  - ➡ The settings are stored.

#### Service note

The menu is used to set whether a service message should be displayed before the service interval expires.

The menu shows the number of days until the next service has to be carried out. If the service is overdue, a message appears.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Service note'* menu is highlighted in blue.
- **<u>2.</u>** Press [*Rotary, push button*].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired setting is selected.
- 4. Press the [BACK] button.
  - ➡ The settings are stored.

#### Visual signal

The menu option is selected so that a visual signal is displayed after centrifugation is complete. The visual signal is the gradual increase in screen brightness.

Opening the lid or pressing any button ends the signal.

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

- **1.** Rotate *[Rotary, push button]* until the *'Visual signal'* menu is high-lighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- 3. Rotate [Rotary, push button] until the desired setting is selected.
- 4. Press the [BACK] button.
  - ➡ The settings are stored.

#### Audible signal

The menu is used to set whether a signal sounds after centrifugation is completed and after warnings, and how loud the signals are.

#### Audible signal after centrifugation is complete

Three consecutive tones sound: "Low", "Medium" and "High".

- The tone sequence is repeated three times.
- This is followed by a pause of three seconds.
- The audible signal is repeated.

The signal is switched off by opening the lid or pressing any button.

#### Audible signal for warnings

Two consecutive tones sound: "Low" and "High".

The audible signal is repeated.

The signal is switched off by opening the lid or pressing any button.

#### Audible signal for errors

The audible signal for errors cannot be configured. If an error occurs, the audible signal sounds at maximum volume.

- A high-pitched sound is emitted.
- The audible signal is repeated.

The signal is switched off by opening the lid or pressing any button.

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Audible signal'* menu is high-lighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired setting is selected at the end of centrifugation.
- 4. Press [Rotary, push button].
  - The next setting is highlighted in blue.
- 5. Rotate [*Rotary, push button*] until the desired setting for warnings is selected.

- 6. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue.
- 7. Rotate [Rotary, push button] until the desired setting for warnings is selected.
- 8. Press the [BACK] button.
  - ➡ The settings are stored.

#### **Display animations**

Display animations are enabled and disabled in the menu.

Display animations are the transitions between different interfaces. If the setting is active, smooth transitions are displayed.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Display animations'* menu is highlighted in blue.
- **2.** Press [Rotary, push button].
  - Menu is displayed.
- 3. Rotate [Rotary, push button] until the desired setting is selected.
- 4. Press the [BACK] button.
  - ➡ The settings are stored.

#### Screen brightness

The screen brightness is set in the menu.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Screen brightness'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- 3. Rotate [Rotary, push button] until the desired setting is selected.
- 4. Press the [BACK] button.
  - ➡ The settings are stored.

#### Energy saving mode

Energy saving mode is set in the menu.

If the setting is active, the display switches off after the entered time if no button is pressed and no messages or signals appear.

#### Set

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Energy saving mode'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- 3. Rotate [Rotary, push button] until 'Yes' is selected.
- 4. Press [Rotary, push button].
  - Option is displayed for setting the time after which energy saving mode is activated.
- 5. Rotate [*Rotary, push button*] until the desired time is set.
  - The time can be set in 30-minute increments up to 8 hours.
- 6. Press the [BACK] button.
  - ➡ The settings are stored.

# Switching off

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Energy saving mode'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- 3. Rotate [Rotary, push button] until 'No' is selected.
- **4.** Press the *[BACK]* button.
  - ➡ The settings are stored.

#### Time

The menu is used to select the time format and set the time.

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Time'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired time format setting is selected.
- 4. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.

- 5. Rotate [Rotary, push button] until the desired value for the hours is selected.
- 6. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.
- 7. Rotate *[Rotary, push button]* until the desired value for the minutes is selected.
- **8.** Press the *[BACK]* button.
  - ➡ The settings are stored.

Date

The menu is used to select the date format and set the date.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Date'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired date format setting is selected.
- 4. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.
- 5. Rotate *[Rotary, push button]* until the desired value for the day or year is selected.
- 6. Press [Rotary, push button].
  - ➡ The next setting is highlighted in blue and displayed in bold.
- 7. Rotate *[Rotary, push button]* until the desired value for the month is selected.
- **8.** Press [*Rotary, push button*].
  - ➡ The next setting is highlighted in blue and displayed in bold.
- **9.** Rotate *[Rotary, push button]* until the desired value for the year or day is selected.
- 10. Press the [BACK] button.
  - The settings are stored.

#### Temperature unit

The "Temperature unit" setting is only available for Mikro 2.0 R.

The menu is used to set whether the temperature is displayed in degrees Celsius or Fahrenheit.

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

- **1.** Rotate *[Rotary, push button]* until the *[Temperature unit]* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired setting is selected.
- **4.** Press [Rotary, push button].
  - ➡ The temperature unit is selected.
- 5. Press the [BACK] button.
  - ➡ The settings are stored.

System language

The menu is used to change the system language.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- 1. Rotate [Rotary, push button] until the 'System language' menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired language is highlighted in blue.
- 4. Press [Rotary, push button].
  - ➡ The language is selected.
- **5.** Press the *[BACK]* button.
  - ➡ The settings are stored.

Changes take effect after a restart.

#### Authorisation level

Some functionalities can only be performed with a sufficiently high authorisation level.

The menu is used to change the authorisation level.

#### For authorisation level 1

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Authorisation level'* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - ➡ Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until Authorisation Level 1 is highlighted in blue.

- 4. Press CONTEXT BUTTON, TOGGLE BUTTON.
  - Context menu is displayed.
- 5. Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
- 6. Press [Rotary, push button].
  - The program is selected.

Any number of programs can be selected.

- 7. Press the [BACK] button.
  - ➡ The settings are stored.
- 8. Press the *BACK* button.
  - ➡ The settings are stored.

#### For authorisation levels 2 and 3

#### Personnel:

- Authorisation level 2
- Authorisation level 3
- 1. Rotate [*Rotary, push button*] until the 'Authorisation level' menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the authorisation level is highlighted in blue.
- **4.** Press [*Rotary, push button*].
  - ➡ The authorisation level is selected.
- 5. Press the [BACK] button.
  - ➡ The settings are stored.

#### Changing the PIN

The PIN for enabling the user levels is changed in the menu.

- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Changing the PIN'* menu is highlighted in blue.
- **2.** Press [Rotary, push button].
  - Menu is displayed.
- **3.** Rotate *[Rotary, push button]* until the 8 digits of the old PIN have been selected.
  - ➡ ✓ symbol is highlighted in blue.
- 4. Press [Rotary, push button].
  - Menu is displayed.
- 5. Rotate [*Rotary, push button*] until the 8 digits of the new PIN have been selected.
  - ➡ ✓ symbol is highlighted in blue.

- 6. Press [Rotary, push button].
  - Menu is displayed.
- 7. Rotate [Rotary, push button] until the 8 digits of the new PIN have been reselected.
  - ➡ ✓ symbol is highlighted in blue.
- 8. Press [Rotary, push button].
  - ➡ The settings are stored.
    - Information line indicates successful execution of the action.

If you forget your PIN, you can reset it using the advanced settings and an authentication file available from Service.

#### Overwriting rotor information

In the menu, information from a previously used rotor is replaced by information from the newly inserted rotor.

The function is available when memory locations for the rotor information are full. This is signalled by warning message 2.2.300.018.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *[Overwriting rotor information]* menu is highlighted in blue.
- 2. Press [Rotary, push button].
  - ➡ The list of rotors is displayed.
- **3.** Rotate *[Rotary, push button]* until the desired rotor is highlighted in blue.
- 4. Press [Rotary, push button].
  - Confirmation query appears.
- 5. Rotate [Rotary, push button] until symbol 🖉 is highlighted in blue.
- 6. Press [Rotary, push button].
  - ➡ Rotor information is overwritten

Information line indicates successful execution of the action.

- 7. Press the [BACK] button.
  - ➡ The submenu is exited.

#### **Default settings**

The menu allows all settings to be reset to their default settings.

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3
- **1.** Rotate *[Rotary, push button]* until the *'Default settings'* menu is high-lighted in blue.

- 2. Press [Rotary, push button].
  - Menu is displayed.
- 3. Press [Rotary, push button].
  - Confirmation query appears.
- **4.** Rotate *[Rotary, push button]* until symbol  $\oslash$  is highlighted in blue.
- 5. Press [Rotary, push button].
  - ➡ User settings are reset to their defaults.
  - ➡ Information line indicates successful execution of the action.
- 6. Press the [BACK] button.
  - ➡ The submenu is exited.

# 6.12 Open source licences

The open source licences used can be displayed in the menu.

#### Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'Open source licences' submenu is selected in the settings menu.

- **1.** Rotate *[Rotary, push button]* to scroll up and down through the open source licences.
- 2. Press the *[BACK]* button.
  - ➡ The submenu is exited.

# 7 Maintenance, cleaning and care

# DANGER

Risk of contamination

Inadequate cleaning or failure to observe the cleaning instructions can lead to contamination risks.

- Observe national and local regulations on safety and accident prevention.
- Observe cleaning instructions.
- Wear personal protective equipment when cleaning the device.
- Comply with laboratory regulations (e.g. TRBAs, the German Protection against Infection Act, hygiene plan) for handling biological agents.

Instructions for cleaning and disinfection

- The device and its accessories must not be cleaned in dishwashers.
- Only perform hand cleaning and liquid disinfection.
- The water temperature must not exceed 25 °C.
- To prevent any corrosion due to use of detergents or disinfectants, it is essential to follow the special application instructions provided by the manufacturers of the detergent or disinfectant.

#### Disinfectant

- Surface disinfectant (not disinfectant for hands or instruments)
- Ethanol as the sole active substance. Do not use an ethanol-propanol mixture to disinfect the viewing window in the lid of the device.
- Concentration is not less than 30%
- pH: 6 8
- Non-corrosive

# 7.1 Cleaning

Maintenance work	Interval in operating hours (h)	To be performed by
Cleaning the device	8	Trained person
Cleaning biosafety systems	100	Trained person
Cleaning accessories	100	Trained person

#### Cleaning the device

- 1. Den the lid.
- 2. Switch off the device and disconnect it from the power supply.
- 3. Remove accessories.
- **4.** Clean the centrifuge housing and the centrifuging chamber with soap or a mild detergent and a damp cloth.
- 5. Remove any detergent residues with a damp cloth after using detergents.
- 6. The surfaces must be dried immediately after cleaning.
- **7.** Dry the centrifuging chamber with an absorbent cloth if condensation forms.

#### Cleaning biosafety systems

- **1.** Clean the biosafety system using the detergent and a damp cloth.
- **<u>2.</u>** Remove any detergent residues with a damp cloth after using detergents.
- 3. Dry the accessories immediately after cleaning using a lint-free cloth and oil-free compressed air. Dry all cavities completely using oil-free compressed air.
- 1. Clean the accessories using the detergent and a damp cloth.
- **2.** Remove any detergent residues with a damp cloth after using detergents.
- 3. Dry the accessories immediately after cleaning using a lint-free cloth and oil-free compressed air. Dry all cavities completely using oil-free compressed air.

# 7.2 Disinfection

**Cleaning accessories** 

Disinfection must always be preceded by cleaning of the components concerned.

Disinfectant concentration and application time according to the manufacturer's instructions.

Maintenance work	Interval in operating hours (h)	To be performed by
Disinfecting the device	if required	Trained person
Disinfecting the accessories	if required	Trained person
Autoclaving	if required	Trained person

CAUTION

### Disinfecting the device

# Risk of electric shock

Liquids that penetrate the device may cause electric shocks.

- The device must be protected from external contact with liquids.
- Do not pour any liquids inside the device.
- The original transport packaging must be used when transporting the device.

### **1.** Open the lid.

- **2.** Switch off the device and disconnect it from the power supply.
- **3.** Remove accessories.
- 4. Clean the housing and centrifuging chamber using disinfectant.
- 5. Remove any disinfectant residues with a damp cloth after using disinfectants.
- **6.** The surfaces must be dried immediately after cleaning.

#### Disinfecting the accessories

- **1.** Disinfect the accessories using the disinfectant.
- **2.** Wet all cavities with bubble-free disinfectant.
- **3.** Remove the disinfectant residues or leave them to dry after using disinfectants.

Autoclaving

The following accessories can be autoclaved at 134  $^{\circ}\text{C}$  / 273  $^{\circ}\text{F}$  2 bar (10 min) up to 10 times:

- Aluminium angle rotors
- Lid

Before autoclaving, the rotor lid must be removed and the seals removed from the rotor.

No statement can be made about the resulting degree of sterility.

Autoclaving accelerates the ageing of materials. It may cause changes to colours.

- **1.** Remove the lid of the rotors before autoclaving.
- 2. After autoclaving, visually check rotors and accessories for damage.
- **3.** Replace damaged parts.

- **4.** Replace the sealing ring if it is cracked, brittle or worn.
- 5. For biosafety systems, replace the sealing rings after autoclaving.

# 7.3 Maintenance

There is a risk of damage to the cooling circuit. The cooling circuit must not be damaged.

Maintenance work	Interval in operating hours (h)	To be performed by
Checking the accessories	100	Qualified mechanics
Checking the biosafety system	100	Qualified mechanics
Centrifuging chamber damage inspection	Annually	Qualified mechanics
Accessories with a limited service life	if required	Qualified mechanics
Replacing centrifuge tubes	if required	Qualified mechanics

WARNING

Checking the accessories

The accessories are to be checked for wear and corrosion damage.
 Remove accessories from operation if they show signs of wear.

- **2.** Check that the rotor is firmly seated.
- Checking the biosafety
   1. Visually check all parts of the biosafety system for damage.

   system
   2. Check the correct installation position of the sealing ring(s) of the biosafety system.
  - 3. Replace the damaged parts of the biosafety system.
  - **4.** Replace the sealing ring if it is cracked, brittle or worn. Replace lids with fixed sealing rings as a complete unit.

Centrifuging chamber damage inspection

Accessories with a limited service life

Do not use accessories if the maximum number of running cycles or the expiry date has been reached.

- The maximum permissible number of running cycles or the expiry date must be complied with.
- The centrifuge is equipped with a cycle counter.

Check the centrifuging chamber for damage.

#### Replacing centrifuge tubes

# CAUTION

Risk of injury

Glass breakage and contaminated liquids in the centrifuge.

- Wear personal protective equipment.
- Stop the centrifuge immediately and disconnect it from the power supply.
- No contact with broken glass or spilled substances.
- Clean the area in accordance with the safety and disposal regulations, watching out for sharp objects and biological material.
- Observe national and local environmental protection and disposal regulations to ensure proper disposal.

Broken parts of the tube, glass splinters and spilled centrifuge material must be removed completely in the event of leakage or if a centrifuge tube breaks. Glass splinters that are not removed will cause further glass breakage.

The rubber inserts and the plastic sleeves of the rotors must be replaced after a glass breakage.

Disinfection must be carried out if the material is infectious.

#### Rotor sealing ring replacement

Fig. 14: Rotor sealing ring positions

- 1 Inner seal
- 2 Outer seal: Lip seal

Remove inner seal and reinstall

# Personnel:

- Trained user
- 1. Remove the inner seal from the groove. Do not use pointed or sharp objects.
- 2. Check the groove for cleanliness and damage.
- 3. Check the seal for damage before inserting it.
- **4.** Insert the seal into the groove.

Remove outer seal (lip seal) and reinstall

- Trained user
- 1. Pull the lip seal out of the groove.
- 2. Check the groove for cleanliness and damage.

- 3. Check the lip seal for damage before inserting it.
- **4.** Align the lips of the lip seal upwards and insert smoothly and without creases into the groove.

# 8 Faults

# 8.1 Fault description

Personnel:

Trained person

In the event of faults that are not listed, contact customer service. Specify the centrifuge type and serial number from the centrifuge rating plate.

Error code number	Cause	Remedy
1.1.001.000 - 1.1.002.004	Error in the tachometer system	Wait until the centrifuge comes to a standstill. Then open the lid. Then switch off the centrifuge. Rotate the rotor by hand and switch on the centrifuge. If the error message cannot be deleted, contact Service.
1.1.004.001 - 1.1.004.005	Lid closure error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.005.000 - 1.1.005.003	Error due to overspeed	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.006.000	Lid closure time-out	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.007.001, 1.1.007.002	Lid closure control error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.010.001 - 1.1.010.005	Error in rotor detection	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.011.001	Power supply interruption	Clear the error by performing a power reset.
1.1.011.002	Mains voltage interruption detected at standstill	Clear the error by performing a power reset. If the fault occurs again, undertake mains voltage stabilising measures. If the mains voltage is stable, contact Service.
1.1.012.001, 1.1.012.002	Invalid machine version	Contact service.
1.1.012.003	Invalid machine component DFC	Contact service.

Error code number	Cause	Remedy	
1.1.012.004	Invalid machine component CFC	Contact service.	
1.1.012.005	Invalid machine component DFC	Contact service.	
1.1.012.006	Invalid machine component CFC	Contact service.	
1.1.014.002	Target speed exceeded for longer than 10 s	Clear the error by performing a power reset. If this occurs repeat- edly, contact Service.	
1.1.014.003	Actual speed too low	Clear the error by performing a power reset. If this occurs repeat- edly, contact Service.	
1.1.025.001	Internal error in the centrifuge con- trol unit	Clear the error by performing a power reset. If this occurs repeat- edly, contact Service.	
1.1.025.002 - 1.1.025.006	Internal error in the centrifuge con- trol unit	Clear the error by performing a power reset. If this occurs repeat- edly, contact Service.	
1.1.026.000 - 1.1.026.050	Sensor defective	Clear the error by performing a power reset. If this occurs repeat- edly, contact Service.	
1.1.027.000, 1.1.027.001	System test Centrifuge control unit error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.	
1.1.028.001 - 1.1.028.007	Internal error in the centrifuge con- trol unit	Clear the error by performing a power reset. If this occurs repeat- edly, contact Service.	
1.1.030.001, 1.1.030.002	Centrifuge control unit error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.	
1.1.031.000 - 1.1.049.003	Internal error in the centrifuge con- trol unit	Clear the error by performing a power reset. If this occurs repeat- edly, contact Service.	
1.1.051.000 - 1.1.055.000	Error in the cooling system	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.	
1.1.057.000	Temperature control failed	The manufacturer recommends compliance with the tempera- ture control specifications. If this occurs repeatedly, contact Service.	
1.1.058.000 - 1.1.058.007	Error in the cooling system	Clear the error by performing a power reset. If this occurs repeat- edly, contact Service.	

Error code number	Cause	Remedy
1.1.060.000	Drive frequency converter error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.060.001	DFC error due to closure	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.061.000, 1.1.061.001	DFC error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.062.000, 1.1.062.001	CFC error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.063.000 - 1.1.063.003	DFC error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.063.004	CFC error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.070.000, 1.1.070.001	Internal error in the centrifuge con- trol unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.071.000, 1.1.071.001	Stop/Open button defective	A defective stop/open button was detected after switching on the centrifuge. Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.086.000 - 1.1.086.012	Internal error in the centrifuge con- trol unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.091.000	Imbalance sensor X-axis defective	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.091.001, 1.1.091.002	Imbalance sensor X-axis offset error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.092.000	Imbalance sensor Y-axis defective	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.092.001, 1.1.092.002	Imbalance sensor Y-axis offset error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.

Error code number	Cause	Remedy
1.1.093.000	Temperature sensor in imbalance sensor defective	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.096.000	Rotor tachometer failure during rotor detection	Delete the error message: Wait until the centrifuge comes to a standstill. Then open the lid. Then switch off the centrifuge. Rotate the rotor by hand and switch on the centrifuge. If the error mes- sage cannot be deleted, contact Service.
1.1.096.001 - 1.1.096.004	Error in the tachometer system	Delete the error message: Wait until the centrifuge comes to a standstill. Then open the lid. Then switch off the centrifuge. Rotate the rotor by hand and switch on the centrifuge. If the error mes- sage cannot be deleted, contact Service.
1.1.097.000, 1.1.098.000	Sensor defective	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.521.001	Communication error, control unit and drive	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.521.003	NAK for communication between control unit and drive	Contact service.
1.1.521.004	No communication between con- trol unit and drive	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.521.005, 1.1.521.006	Communication error drive control unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.400.001	Stop/Open button malfunction detected	Contact service.
2.1.420.001	Centrifuge control unit type error	Switch off the device. Contact service.
2.1.501.001	Centrifuge control unit illogical information	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.541.001	Power supply interruption	Clear the error by performing a power reset.
2.1.551.001	Error message logging	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.

Error code number	Cause	Remedy
2.1.600.001	Error in the application	This error message can only be deleted by performing a power reset.
2.1.620.000	ServiceAuth file incorrect	Clear the error by performing a power reset.
2.1.800.002 - 2.1.800.010	Error message logging	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.901.002	Drive is not ready	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.901.003	Function test not possible	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.901.004	Centrifuge control unit version	Contact service.

2.1.901.004	Centrifuge control unit version error	Contact service.
2.1.901.005	Centrifuge control unit configura- tion error	Contact service.

# 8.2 Warning descriptions

Personnel:

Trained person

In the event of faults that are not listed, contact customer service. Specify the centrifuge type and serial number from the centrifuge rating plate.

Warning code number	Cause	Remedy
1.2.003.001	The centrifuge has detected an imbalance.	Please open the lid when the cen- trifuge is at a standstill. Check loading. Ensure even weight distribution.
1.2.011.000	A power failure occurred during centrifugation.	Open the lid when the centrifuge is at a standstill. The manufacturer recommends checking the result of the centrifugation.
2.2.300.001	The centrifuge has detected a rotor change.	Acknowledgement of the warning only possible at standstill. The centrifuge will check or adjust the current centrifugation parame- ters for the current rotor.

Faults

Warning code number	Cause	Remedy
2.2.300.002	Starting of centrifugation without a rotor.	Acknowledgement of the warning only possible at standstill. The centrifuge has detected that centrifugation has been started without a rotor. Only start centrifu- gation with a rotor.
2.2.300.003	The permissible number of cycles for the rotor currently in use has been reached.	The rotor currently in use must be replaced immediately with a new rotor. We warn against further use. Further use only at your own risk.
2.2.300.004	The permissible number of cycles for the rotor currently in use has been reached.	Use of the current rotor poses a safety risk. Use only at your own risk.
2.2.300.005	The centrifuge has detected a rotor change.	The centrifuge will check or adjust the current centrifugation parame- ters for the current rotor.
2.2.300.010, 2.2.300.011	The permitted number of centri- fugations has been reached in spite of a faulty rotor identification system.	Further centrifugation is not pos- sible. Contact service.
2.2.300.012	The rotor identification system (tag) is faulty or no rotor is fitted.	Only a limited number of centrifu- gations are possible in this state. It is strongly recommended that you contact Service if the rotor identification system malfunctions. If no rotor is fitted, the start of centrifugation is terminated with a warning message.
2.2.300.013	The permissible operating period for the rotor has been exceeded.	The rotor currently in use must be replaced immediately with a new rotor. We warn against further use. Further use only at your own risk.
2.2.300.014	The rotor identification system (reader) is faulty.	Only a limited number of centrifu- gations are possible in this state. Contact service.
2.2.300.015	The rotor identification system reports an error.	A limited number of centrifugations are still possible. Contact service. More information on the number of centrifugations at the start of this.
2.2.300.016	The current rotor is not approved for this centrifuge. It is not possible to start centrifugation.	Approved rotors can be found in the operating instructions. Contact service.

**Ilettich** 

Warning code number	Cause	Remedy
2.2.300.017	The current rotor does not comply with the Hettich standard.	It is not possible to start centrifu- gation. Please remove this rotor from the centrifuge immediately. Contact service.
2.2.300.018	The memory locations for rotor information are full. It is currently not possible to start centrifugation.	Select an entry in "Overwrite rotor info" that may be overwritten with the data of the current rotor. This action is only possible when the new rotor is in the centrifuge.
2.2.300.019, 2.2.300.020	The permitted number of centri- fugations has been reached in spite of a faulty rotor identification system.	The manufacturer recommends reactivating the centrifugation allo- cation. Contact service.
2.2.310.001, 2.2.310.002	The user has cancelled the cur- rent centrifugation by pressing the Stop button.	Acknowledgement of the warning only possible at standstill. The manufacturer recommends checking the result of the centrifu- gation.
2.2.320.001	User has cancelled the data export.	Remove USB flash drive or restart data export.
2.2.340.001	The desired temperature is above the current temperature. This cen- trifuge can only cool.	To achieve a higher temperature than the current one, carry out a standard centrifugation.
2.2.550.013	The update package does not contain a more up-to-date soft-ware version for the centrifuge.	The software update is cancelled.
2.2.550.014	The software version of the HBI update file was not accepted.	Check the version of the software update used.
2.2.550.015	The ZS software version has not been updated.	Repeat software update.
2.2.560.001	Failed to mount the USB flash drive.	Use another USB flash drive. If this occurs repeatedly, contact Service.
2.2.560.002	Unmounting the USB flash drive failed.	Remove the USB flash drive and restart the centrifuge. If this occurs repeatedly, contact Service.
2.2.560.003	The system reports: "No USB flash drive available".	Remove the USB flash drive and insert it again. Use a different USB flash drive if the warning message appears again. If there is a per- manent warning message, contact Service.

Warning code number	Cause	Remedy	
2.2.560.004	Insufficient storage space on USB flash drive for data export.	Use a USB flash drive with suffi- cient free storage space. Recom- mendation: at least 8 GB	
2.2.560.005	Failed to write to USB flash drive. Data may not have been com- pletely written to the USB flash drive.	Check USB flash drive and restart export.	
2.2.560.006	The file system of the USB flash drive does not correspond to the system specifications. Data storage is only possible with the file systems: FAT, FAT32, ext2, ext3, ext4.	Use a USB flash drive with a supported file system.	
2.2.560.007	Deletion of a file on the USB flash drive has failed. Copying process was cancelled.	Clean up the USB flash drive and export again.	
2.2.560.008	Remounting the USB flash drive has failed. Files may not have been copied completely.	Check USB flash drive and per- form the export again.	
2.2.570.001	Next service imminent.	Contact service.	
2.2.600.002	A defective database was detected and restored.	Acknowledge warning message. If this occurs repeatedly, contact Service.	
2.2.600.003	The database restore has failed. The database has been recreated. Programs and settings are no longer available.	Contact service.	
2.2.600.004	Error when deleting the system logs.	Perform "Reset to delivery status" again. If this occurs repeatedly, contact Service.	
2.2.600.010	A problem with the date or time was detected when the centrifuge was started.	Please correct the date or time in centrifuge. If this occurs repeatedly, contact Service.	
2.2.600.011	The battery of the internal clock is discharged. The service request was activated out of schedule.	To replace the internal battery, contact Service.	
2.2.600.021	Machine version rejected.	Limited functionality, no centrifuga- tion possible. Contact service.	
2.2.600.040	Balancing of the motor and DFC is mandatory.	Centrifugation is only possible after balancing. Contact service.	

# 8.3 Performing a MAINS RESET

- **1.** Set the mains switch to [0].
- 2. Wait 10 seconds.
- 3. Set the mains switch to [/].

## 8.4 Emergency release

The lid cannot be unlocked by the motor in the event of a power failure. Emergency unlocking by hand must be performed.

#### 

#### Danger to life due to electric current

Contact with live parts can be fatal.

- Work on electrical devices must only be carried out by a qualified electrician.
- Work must only be carried out on disconnected devices.
- Check that disconnected devices are de-energised before carrying out any work.
- Some device components may not be de-energised even after being switched off. These are specially marked. Work on these components must only be carried out according to the instructions on the label.
- Only insulated tools may be used when working on the electrical system.
- Observe national and local regulations on safety and accident prevention.

#### WARNING

#### Risk of cuts and crushing

Risk of injury due to moving rotor.

 Do not open the lid until the rotor has come to a complete stop.

#### Personnel:

- Trained user
- **1.** Look through the window in the lid to ensure that the rotor is stationary.
- **2.** Insert an Allen key horizontally into the hole (1).
- **3.** Rotate the Allen key anticlockwise until the lid opens.

1 Hole

# 8.5 Replacing the mains input fuse

## 

#### Danger to life due to electric current

Contact with live parts can be fatal.

- Work on electrical devices must only be carried out by a qualified electrician.
- Work must only be carried out on disconnected devices.
- Check that disconnected devices are de-energised before carrying out any work.
- Some device components may not be de-energised even after being switched off. These are specially marked. Work on these components must only be carried out according to the instructions on the label.
- Only insulated tools may be used when working on the electrical system.
- Observe national and local regulations on safety and accident prevention.

#### Personnel:

- Trained user
- The mains fuses are located next to the mains switch.

The mains switch is in switch position [O]

- **1.** Disconnect the mains cable from the device plug.
- **2.** Press the snap locks (2) against the fuse holder (1) and pull them out.
- **3.** Replace the defective mains input fuses.

Only use fuses with the nominal value specified for the type: see the table below.

- **4.** Push in the fuse holder (1) until the snap lock engages.
- 5. Reconnect the device to the mains.

Model	Туре	Fuse	Order no.
MIKRO 2.0	2404 2407	T8 AH/250V	E738
MIKRO 2.0 R	2406 2408	T8 AH/250V	E738

# 9 Disassembly

Disassembly is carried out in reverse order to assembly.

#### 10 Spare parts

Only use original spare parts from the manufacturer, approved accessories and wear parts.

Fuse holder
 Snap lock

# 11 Storage

- Insert property damage reference
   The minimum room floor area for storage must correspond to the same minimum room floor area for operation.
- Do not damage the base body and protruding components when setting the device down.
- Store the device in its original packaging.
- Use a wooden, rubber or plastic base.
- Clean thoroughly and preserve before prolonged storage.
- Protect against dust and coarse dirt.
- Store in a closed, dry room.
- Observe storage and ambient conditions.

Specification	Value
Permissible temperature range	-20 to 60 °C
Relative humidity (non-condensing)	10 to 80%

# 12 Disposal

#### 12.1 General instructions

<ul> <li>The device can be disposed of via the manufacturer.</li> <li>A Return Material Authorisation (RMA) form must always be requested for a return.</li> <li>If necessary, contact the Technical Service Department of the manufacturer:</li> <li>Andreas Hettich GmbH</li> <li>Föhrenstraße 12</li> <li>78532 Tuttlingen, Germany</li> <li>Phone: +49 7461 705 1400</li> </ul>
- E-Mail: service@hettichlab.com
The cooling circuit must only be dismantled by qualified personnel.
WARNING
Risk of soiling and contamination
 Damage to the environment and health due to improper disposal.
<ul> <li>Observe national and local environmental protection and disposal regulations to ensure proper disposal or recy- cling. Separate metals, non-metals, composite materials</li> </ul>

- and auxiliary materials by type and dispose of them in an eco-friendly manner.
- The device may only be disassembled and disposed of by trained and authorised service technicians.

The device is intended for the commercial sector ("Business to Business" - B2B).

According to Directive 2012/19/EU, the devices may no longer be disposed of with household waste.

The devices are assigned to the following groups according to the Stiftung Elektro-Altgeräte Register (EAR (German foundation under civil law)):

Group 5 (small devices)

The crossed-out wheelie bin symbol indicates that the device must not be disposed of with household waste. Regulations governing disposal of such devices may differ in individual countries. If necessary, contact the supplier.

13 Glossary	
Autoclaving	Autoclaving is a process for sterilising materials and objects using pressur- ised saturated steam in an autoclave. An autoclave is a closed device capable of generating high pressure and high temperatures to kill microor- ganisms such as bacteria, viruses, fungi and spores.
Basic-UDI-DI	The basic version of the Unique Device Identification, which provides a lifelong identification feature for medical devices within the EU.
Brake level	Determines how quickly the centrifuge is decelerated from maximum speed to a complete standstill, influenced by various levels from gentle to fast.
Centrifugation	The process of applying centrifugal force to separate components of a solution based on their density.
Centrifugation allocation	Allocation (number) of approved centrifugations that the customer may per- form despite a malfunction in the rotor identification system.
Centrifuge lid	The part of the centrifuge that remains closed during operation to ensure safety and protect the integrity of the samples.
Commissioning	The process of setting up and starting the centrifuge before first use or after maintenance.
G-force (RCF – Relative Centri- fugal Force)	The force exerted on a sample in the centrifuge, expressed as a multiple of the acceleration due to gravity.
General information about the glossary	* This glossary provides a comprehensive overview of key terms that are important when using the centrifuge and helps laboratory personnel to better understand the documentation and operating instructions for the centrifuge.
PIN	Personal identification number used to access secure centrifuge functions.
Power reset	A procedure for resetting the electrical supply to the centrifuge in order to correct system errors or restart the system.
Ramp-up level	Defines the rate of acceleration of the centrifuge up to the target speed. Different levels enable gentle or rapid acceleration depending on sample sensitivity.
RCF (Relative Centrifugal Force)	The force acting on the sample during centrifugation, measured as a mul- tiple of the acceleration due to gravity (g).
Residual current circuit breaker (RCD)	The residual current circuit breaker interrupts the power supply in the event of a residual current and thus protects against electric shocks and fires.
Risk of contamination	Risk of contamination of the samples or the working environment, in particular by pathogenic organisms or chemical substances.
Rotor	The rotating part of the centrifuge in which the sample containers are inserted. Rotors can vary depending on the application.
Safety area	The area around the centrifuge that must be free of obstacles to ensure safety during operation.
UDI (Unique Device Identifica- tion)	A worldwide system for the unique identification of medical devices over their entire service life.

# Tlettich

#### Index 14

# Α

Accessories	
autoclaving	60
biosafety systems	59
checking	61
disinfection	60
Authorisation level	55
Authorisation levels	
Authorisations	7
Autoclaving	60
_	

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Biosafety	system
-----------	--------

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Centrifugation	
Short-term centrifugation	37
with higher substance density	28
with time preselection	36
Centrifuge tubes	
filling	26
replacing	62
Certifications.	13
Changing the PIN	56
Cleaning	
biosafety systems.	59
Connecting the centrifuge	20
Continuous operation.	35
-	

# D

Date format	54
biosafety systems	59 60
Device information	
Data export to USB flash drive	47
Error history	46
Querying.	46
Rotor info	45
Service contact details	47
System information	45
Display animations.	52
Disposal	73
E	
Energy saving mode	52
F	
Filling	26
G	
General safety instructions	9
L	
Labels	
on the device	15
on the packaging	14

Lid
Automatic start after lid closure.49Lid opens after centrifugation.49Lid opens when switched on.49the lid.23, 24Loading.26Logos.13
Μ
MAINS RESET
menu
Menu
changing.33Main menu.33Program menu.34Quick Adjust menu.34, 39selecting.33Selection of submenus.35Settings menu.35
0
Open source licences
a program link
Operator obligations
Obligations of the operator
Р
Program
changing
retrieving
Program favourite creation
R
Ramp-up level.29Rating plate.12Relative centrifugal force

RCF	27
Removing	
Symbols	26
the rotor	27
Return.	15
Rotor	
installation	24
removal	24
Rotor information.	57
Rotor lid	
fitting	25
removal	25
Runtime	29
changing 40, 40, 40, 40, 40, 40, 40, 40, 40, 40,	42
Runtime display	48

# s

Safety instructions	9 17 52 50 19
Signal         Audible signal.         Visual signal.         Speed.       39,         Starting point in Quick Adjust.         Step size for runtime settings.         Storage conditions.         Switching off.         Switching off.         Switching the unit on.         Switching on for the first time.         Symbols.       5,         System language.         T	51 50 42 47 48 23 22 21 31 55
Temperature unit	54 53 17
removal.	18
Unpacking	19

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Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)

 (a) Observe the tube manufacturer's instructions.
(a) nonformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance" for bio safety systems in chapters indies on sa and servicing". 5) recommended for high-speed centrifugation 6) Load only each second position of the rotor 7) RZB / RCF top row 8) RZB / RCF bottom row